



MAGNETIC STIRRERS



2mag MAGNETIC STIRRERS WELL-PROVEN CONCEPTS - INNOVATIVE PRODUCTS



WELL-PROVEN CONCEPTS - INNOVATIVE PRODUCTS



- 100% wear- and maintenance-free
- 30% more stirring power than comparable systems
- Super-flat, space-saving stirring drives ideal for robotics
- Made for permanent operations
- Heat resistant and submersible
- Synchronous speed at all stirring positions
- Zero-stand-by power supply units
- Individual configurations fast and economic realizable
- Experience with inductive and magnetic drive concepts since 1974

magnetic ^emotion



INDUCTIVE DRIVE -MAINTENANCE- AND WEAR-FREE

Our magnetic drives are working on the inductive principle. This means that instead of conventionally problematic motors with pull magnets, 2mag is working with stationary coils, which are defined flown through by alternating currents. The thereby generated electromagnetic rotating field is 100% wear- and maintenance-free.

The drive does not have any mechanically moved parts, as e.g. belts, bearings or motors. Therefore, the drives are reliable and absolutely wear-free even at high operation temperatures (up to +300 °C).

2mag MAGNETIC STIRRERS WELL-PROVEN CONCEPTS - INNOVATIVE PRODUCTS



Single stirring point Multiple stirring points Up to 3 litres stirring volume, without motor MX 1 eco accuMX accuMX IwMX accuMX WMXdrive 1 XS (HT) MX 15 eco WX 1 for 0 10 litres stirring volume, without motor MX 1 S MX 1 MX 4 MS MX 4 MS MX 6 MX 1 XL MX 4 MS MX 4 MX 6 MX 1 S MX 1 XL MX 4 MS MX 4 MX 6 MX 1 S MX 1 XL MX 4 MS MX 1 XL MX 4 MS MX 1 XL MX 4 MS MX MY 1 XL MX MX 1 MU MX MX 1 MI MX MX 1 MU			
MX 1 eco accuMX LUMX CuvetteMXdrive 1 MXdrive 1 XS (HT) MXdrive 1 eco (HT) MX 15 eco Up to 10 litres stirring volume, without motor MX 1 MX 1 MX 4 MS MX 4 MS MX 4 MS MX 6 MX 15 MX 6 MX 15 MX 8 XL MX 15 MX 12 XL MX drive 1 (HT) atexMX drive MXdrive 6 (HT) MX drive 15 (HT) MX drive 15 (HT) MX drive 6 (HT) MX drive 6 (HT) MX drive 6 (HT) MX drive 15 (HT) MX drive 15 (HT) MX drive 15 (HT) MX drive 15 (HT) MX drive 10 For cell cultures, warming-free bioMIX drive MX drive 1 MX 1 MX drive		Single stirring point	Multiple stirring points
accuMIX IUM		Up to 3 litres s	tirring volume, without motor
IUMIX w cuvetteMIXdrive 1 w MIXdrive 1 XS (HT) w MIXdrive 1 eco (HT) w IUp to 10 litres stirring volume, without motor MIX 1 MIX 4 MIX 6 MIX 6 MIX 8 XL MIX 15 MIX 8 XL MIXdrive 1 (HT) w atexMIXdrive w MIXdrive 6 (HT) w MIXdrive 15 (HT) w MIXdrive 6 (HT) w MIXdrive 6 (HT) w MIXdrive 6 (HT) w MIXdrive 15 (HT) w MIXdrive 6 (HT) w MIXdrive 10 (UT) w MIXdrive 10 (UT) <td></td> <td>MIX 1 eco</td> <td></td>		MIX 1 eco	
CuvetteMIXdrive 1 XS (HT) MIXdrive 1 XS (HT) MIX drive 1 XS (HT) MIX 15 eco Up to 10 litres stirring volume, without motor MIX 1 Seco MIX 1 MS MIX 4 MS MIX 1 S M		accuMIX	
MXdrive 1 XS (HT) MXdrive 1 eco (HT) MX 15 eco Up to 10 litres stirring volume, without motor MX 1 MX 1 MX 4 MS MX 6 MX 6 MX 15 MX 8 XL MX 6 (HT) MX drive 1 (HT) MX drive 1 (HT) MX drive 6 (HT) MX drive 6 (HT) MX drive 6 (HT) MX drive 6 (HT) FABdrive FABdrive For cell cultures, warming-free bioMIX 1 bioMIX 1 bioMIX 1 MX drive 1 (HT) MX drive 1 XL MX drive 1 XL MX drive 2 / 3 / 4 For cell cultures, warming-free bioMIX 1 MX drive 2 / 3 / 4 For microtiter plates MX drive 6 / 12 / 24 / 96 MTP Heatable stirring systems		IuMIX	
MIX drive 1 eco (HT) MIX 15 eco Up to 10 litres stirring volume, without motor MIX 1 MIX 1 MIX 4 MS MIX 6 MIX 6 MIX 8 MS MIX 15 MIX 15 MIX 16 MIX 175 MIX 175 MIX 100 MIX 1010 MIX 112 XL MIX 10100 MIX 10100 MIX 11000 MIX 110000 MIX 110000 MIX 110000 MIX 1100000 MIX 1100000 MIX 1100000000000000000000000000000000000	や	cuvetteMIXdrive 1	
MIX 15 eco Up to 10 litres stirring volume, without motor MIX 1 MIX 1 MIX 4 MS MIX 4 MS MIX 4 MS MIX 4 MS MIX 5 MIX 15 MIX 15 MIX 12 XL MIX 1	や	MIXdrive 1 XS (HT)	
WIX 1 MIX 4 MS MIX 4 MS MIX 6 MIX 6 MIX 6 MIX 10 MIX 6 MIX 11 MIX 6 MIX 2000 MIX 8000 MIX 2000 MIX 12 XL MIX 11/2 MIX 15 (HT) MIX 11/2 MIX 10/2 MIX 11/2 MIX 10/2 MIX 11/2 MIX 10/2 MIX 11/2 MIX 11/2	や	MIXdrive 1 eco (HT)	
MIX 1 MIX 4 MS MIX 4 MS MIX 4 MS MIX 6 MIX 10 MIX 10 MIX 10 MIX 10 MIX 10 MIX 112 XL MIX 12 XL MIX			MIX 15 eco
MX 4 MS MX 6 MX 10 MX 15 MX 8 XL MX 12 XL MX 14rive 6 (HT) MX 14rive 6 (HT) MX 14rive 6 (HT) MX 14rive 7 MX 14rive 8 MX 14rive 9 MX 14rive 9 MX 14rive 9 MX 14rive 10 MX 14rive MX 11 MX 1 MX 1<		Up to 10 litres s	stirring volume, without motor
MIX 6 MIX 15 MIX 8 XL MIX 8 XL MIX 12 X		MIX 1	
MiX 15 MiX 8 XL MiX 12 XL MiX 13 VH MiX 15 (HT) MiX 15 (HT) MiX 15 (HT) MiX 16 (HT) MiX 17			MIX 4 MS
MIX 8 XL MIX 12 XL MIX 12 XL MIX 12 XL MIX 12 XL MIX 10 XC MIX 10 XC MIX 10 XC MIX 10 XC MIX 1 XL MIX 1 X			MIX 6
MX 12 XL MX drive 1 (HT) a texMX drive texMX drive A MX drive 6 (HT) A MX drive 6 (HT) A MX drive 6 (HT) A MX drive 15 (HT) A MX drive 60			MIX 15
MIXdrive 1 (HT) atexMIXdrive atexMIXdrive atexMIXdrive atexMIXdrive atexMIX atexMIX atexMIX atexMIX atexMIX atexMIX aterMIX atexMIX aterMIX			MIX 8 XL
A atexMIXdrive A atexMIXdrive A atexMIXdrive A mixdrive 6 (HT) A mixdrive 15 (HT) A mixdrive 60 (HT) A mixdrive 60 (HT) A mixdrive 1 XL maxMIX Mixdrive 1 XL maxMIX A mixdrive FABdrive A mixdrive For cell cultures, warming-free bioMIX 1 B bioMIX 1 Mixdrive 1 B bioMIX 1 B bi			MIX 12 XL
MIXdrive 6 (HT) MIXdrive 15 (HT) MIXdrive 60 (HT	や	MIXdrive 1 (HT)	
NIX drive 15 (HT) NIX drive 60 (HT) Up to 1,000 litres stirring volume, strong MIX 1 XL NIX 1 XL MIX drive 1 XL maxMIX MAXdrive Katrive FABdrive SteriMIXdrive Image: SteriMIXdrive Image: SteriMIXdrive 1 Image: SteriMIXdrive 2/3/4 Image: SteriMIXdrive 5/12/24/96 MTP Image: SteriMIXdrive 5/12/24/96 MTP	や	atexMIXdrive	
Image: Nit X drive 6 / 12 / 24 / 96 MTP Image: Nit X drive 1 XL Image: Nit X drive Image: Nit X drive 1 Image: Nit X drive 2 / 3 / 4 Image: Nit X drive 6 / 12 / 24 / 96 MTP Image: Nit X drive 6 / 12 / 24 / 96 MTP Image: Nit X drive 6 / 12 / 24 / 96 MTP			ペ MIXdrive 6 (HT)
Image: With X to X			ℵ MIXdrive 15 (HT)
MIX 1 XL R MIX drive 1 XL maxMIX MAXdrive MAXdrive FABdrive FABdrive For cell cultures, warming-free bioMIX 1 bioMIX 1 bioMIX 1 bioMIX 4rive 1 For microtiter plates MIXdrive 6 / 12 / 24 / 96 MTP Heatable stirring systems			≈ MIXdrive 60 (HT)
MIXdrive 1 XL maxMIX MAXdrive MAXdrive MAXdrive FABdrive steriMIXdrive steriMIXdrive maxmanuscup ma		Up to 1,000 l	itres stirring volume, strong
maxMIX MAXdrive MAXdrive FABdrive isteriMIXdrive isteriMIXdrive MiXdrive 1 isteriMIXdrive 1 isteriMIXdrive 1 MiXdrive 2/3/4 isteriMIXdrive 2/3/4 MIXdrive 6/12/24/96 MTP Heatable stirrIng systems		MIX 1 XL	
MAXdrive MAXdrive MAXdrive FABdrive SteriMIXdrive MIXdrive MIXdrive 1	や	MIXdrive 1 XL	
PAB drive Paber in MIX drive Paber in MIX drive Paber in MIX 1 Paber in MIX drive 1 Paber in MIX drive 2 / 3 / 4 Paber in MIX drive 6 / 12 / 24 / 96 MTP Heatable stirring systems		maxMIX	
steriMIXdrive For cell cultures, warming-free bioMIX 1 bioMIX 1 bioMIXdrive 1 bioMIXdrive 2 / 3 / 4 For microtiter plates NIXdrive 6 / 12 / 24 / 96 MTP Heatable stirring systems	や	MAXdrive	
For cell cultures, warming-free bioMIX 1 bioMIXdrive 1 bioMIXdrive 2/3/4 For microtiter plates MIXdrive 6/12/24/96 MTP Heatable stirring systems	や	FABdrive	
bioMIX 1	や	steriMIXdrive	
bioMIXdrive 1 bioMIXdrive 2 / 3 / 4 For microtiter plates MIXdrive 6 / 12 / 24 / 96 MTP Heatable stirring systems		For cell	cultures, warming-free
 bioMIXdrive 2 / 3 / 4 For microtiter plates MIXdrive 6 / 12 / 24 / 96 MTP Heatable stirring systems 		bioMIX 1	
For microtiter plates	や	bioMIXdrive 1	
№ MIXdrive 6 / 12 / 24 / 96 MTP Heatable stirring systems			Normal Karive 2 / 3 / 4
Heatable stirring systems		For	microtiter plates
			浸 MIXdrive 6 / 12 / 24 / 96 MTP
		Heata	ible stirring systems
hotMIX 1		hotMIX 1	
ル STIRRING DRYBATH 15-100			≈ STIRRING DRYBATH 15-100
ル STIRRING DRYBATH 15-250			⋧ STIRRING DRYBATH 15-250
ル STIRRING DRYBATH 8-250 ERL			≈ STIRRING DRYBATH 8-250 ERL
₹ STIRRING HOTPLATE 6			≈ STIRRING HOTPLATE 6
ル STIRRING HOTPLATE 15			≈ STIRRING HOTPLATE 15
with external control unit		a with	external control unit

magnetic ^emotion

The main importance of our passion lies upon our customers and our products! 2mag is providing you more than 30 years of competence in the conception, development, production and sales of magnetic and inductive drive concepts, as e.g. magnetic stirrers. This experience paired with a great passion and emotion for our customers – and our products – enables the fast and economic creation of solutions precisely for your needs.

"Customizing", the manufacturing of special made products, is our business, our competence and our passion! Individual requests will result in individual – and even more important – also economically realizable solutions.

Ask us – test us – we are there for you!

emotion for motion

- Customer-oriented
- Short communication ways competent team
- Direct access to experienced contact persons
- Fast reactions
- Innovative products
- Well-proven concepts
- Experience with inductive and magnetic drive concepts since 1974

TABLE OF CONTENTS



Page

28

29

30 31

32

33

34 35

36 37

38 39

40 41 42

43 45

46

47

48

51

52 54

55

56

magnetic ^emotion



What is the effect of a magnetic North pole without a South pole? What is the effect of an electric + Pole without a - Pole? What is the effect of Techniques without Competence? What is the effect of a Product without Passion? What is the effect of a Supplier without a Customer?

According to experience, always 2 are needed to achieve a successful dialogue and interaction. The main focus of our action and business philosophy are our customers and their needs and requirements.

Δ

Single point stirrers	Page	5.	Cell culture stirrers
with internal control unit MIX 1	5		with internal control unit bioMIX 1
MIX 1 eco accuMIX IuMIX	6 7 8		with external control unit bioMIXdrive 1 / 2 / 3 / 4
MIXphilosophy	9	6.	Control units for bioMIXdrive bioMIXcontrol / bioMIXcontrol S
with external control unit	10		bioMIXcontrol 4MS
MIXdrive 1 MIXdrive 1 eco MIXdrive 1 XS	10 11 12	7.	Microtiter plate stirrers MIXdrive 6 / 12 / 24 / 96 MTP
cuvetteMIXdrive 1	13	8.	Control unit for MIXdrive MTP
Multiple stirrers			
with internal control unit		9.	Industrial stirrers
MIX 6 / MIX 15 MIX 15 eco	14 15		with internal control unit
MIX 13 ECO MIX 8 XL / MIX 12 XL	16		MIX 1 XL maxMIX
MIX 4 MS	17		
MIXphilosophy	18		with external control unit MIXdrive 1 XL / MIXcontrol XL MAXdrive / MAXcontrol
with external control unit MIXdrive 6 / 15 / 60	19		FABdrive / FABcontrol
Control units for MIXdrive			mixWATCH
MIXcontrol eco	20	10.	Specialised stirrers
MIXcontrol eco DINrail	21		steriMIXdrive / steriMIXcontrol
MIXcontrol 20	22		atexMIXdrive
MIXcontrol 40	23		atexMIXcontrol / cleanroom
Heatable stirrers		11.	Accessories
Single point stirrers			Stirring bars
hotMIX 1	24		Stirring bar remover Extension cords / Distributors
Multiple stirrers			Accessories for multiple stirrers Accessories for heatable stirrers
Hotplates STIRRING HOTPLATE 6 / 15	25		Accessories for industrial stirrers Water baths
Drybaths STIRRING DRYBATH 15-100 / 15-250 STIRRING DRYBATH 8-250 ERL) 26 27	13.	Individual stirrers Sustainability Contact

1.

2.

3.

4.

SINGLE POINT STIRRERS WITH INTERNAL CONTROL

Designed for chemical, biotechnical and medical applications of all kinds, developed for the daily lab use and continuous operation.



MIX 1

Ultra flat, wear-free magnetic stirrer with 1 stirring position, up to 10,000 ml stirring volume, also suitable for viscous media.

- *№* Maintenance- and wear-free
- ℵ Up to 10 litres or for viscous media
- ∂ Outstanding long-life quality
- ֎ Developed and made in Germany

magnetic ^emotion



Туре	MIX 1		
Stirring points	1		
Stirring volume/stirring point	1 - 10,000 ml		
Speed range	100 - 2,000 rpm		
Stirring power (max.)	20 W		
Power setting	5/10/15/20 W (4-step)		
Material housing	stainless steel		
Permitted operation conditions	-10 °C up to +50 °C (at 80% humidity)		
Electrical data	100-240 V / 50-60 Hz / 1.5 A		
Dimensions (WxDxH)	188 x 238 x 38 mm		
Protection category	IP65		
Weight (gross)	approx. 4.2 kg		
Order no.	30001		

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

100% maintenance-free and wear-free inductive 2mag-Magnetic-Drive concept for powerful mixing of solutions, speed range from 100 up to 2,000 rpm, jerk-free stirring even at low speeds.

4-step power setting for high power with regard to large and viscous stirring volumes and reduced power for the warming-free continuous operation. Clear digital display for speed- and power settings, SoftStart for reliable catching/centering and safe accelerating of the stirring bar. QuickSet for quick setting of start- and maximum speed, automatic memory function to save the last settings, outstanding long-life cycle of the hermetically sealed drive system, resistant stainless steel housing for easy and fast cleaning.

Accessories

ASTEROID 40 - Stirring bar Stirring bar remover PTFE

SINGLE POINT STIRRERS WITH INTERNAL CONTROL

Designed for chemical, biotechnical and medical applications of all kinds, developed for the daily lab use and continuous operation.



MIX 1 eco

Ultra flat, wear-free magnetic stirrer with 1 stirring position, up to 3,000 ml stirring volume, also suitable for the operation with aggressive media.

∂ Price-optimized

- ∛ Up to 3,000 ml
- ✤ Developed and made in Germany

magnetic ^emotion

1



3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Price-optimized magnetic stirrer, 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept for mixing of solutions, speed range from 120 up to 1,400 rpm, jerk-free stirring even at low speeds.

SoftStart for reliable catching/centring and safe accelerating of the stirring bar, fully encapsulated stainless steel housing, extremely robust construction.

Accessories

ASTEROID 25 - Stirring bar Stirring bar remover PTFE

SINGLE POINT STIRRER FOR MOBILE OPERATION

Designed for the mobile and mains-independent operation, ideal for chemical, biotechnical and medical applications, en route and far away from laboratory and infrastructure.



accuMIX

7

Ultra flat, wear-free magnetic stirrer with 1 stirring position, up to 3,000 ml stirring volume, robust and firm construction for rechargeable battery- and mains operation.

- ∂ Battery operation
- 🗞 Mobile use
- Up to 3,000 ml
- ℵ Robust and durable
- Neveloped and made in Germany





Туре	accuMIX		
Stirring points	1		
Stirring volume/stirring point	1 - 3,000 ml		
Speed range	100 - 1,600 rpm		
Stirring power (max.)	10 W		
Power setting	2/10 W (2-step)		
Material housing	stainless steel		
Permitted operation conditions	-10 °C up to +50 °C (at 80% humidity)		
Electrical data	100-240 V / 50-60 Hz / 1.7 A		
Dimensions (WxDxH)	215 x 130 x 40 mm		
Protection category	IP30		
Weight (gross)	approx. 2.6 kg		
Order no.	30401		

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Mobile magnetic stirrer, 100% maintenance- and wear-free by inductive 2mag-Magnetic-Drive concept for mixing of solutions, speed range from 100 up to 1,600 rpm, jerk-free stirring also at low speeds, two-step power setting for high power with regard to large and viscous stirring amounts and reduced power for battery-saving long-term operation. SoftStart for reliable catching/centering and safe accelerating of stirring bar, stainless steel housing, very robust construction.

Constructed for the independent battery- and stationary mains operation, equipped with 8 commercially available high performance rechargeable batteries (Mignon/AA), universal charging/power supply device, intelligent battery charging station with single controlled charging slots as well as battery service function and 12 V adapter cable for the operation at car sockets respectively cigarette lighter.

Accessories

ASTEROID 40 - Stirring bar Stirring bar remover PTFE

SINGLE POINT STIRRER WITH ILLUMINATED STIRRING PLATE

Designed for chemical, biotechnical and medical applications of all kinds, developed for the hard daily lab use and uncompromising continuous operation. Particularly glare-free illumination of samples for the optical test range, e.g. color changes, titrations ...



IuMIX

Ultra-flat, wear-free magnetic stirrer with 1 large stirring position and a luminescent plate, up to 3,000 mL stirring volume, also suitable for viscous media. Stirring function and illumination function separately switchable.



- ℵ Large stirring position, luminescent
- ℰ Homogenous brightness, variable
- ℵ Maintenance- and wear-free
- Note: Note:
- ֎ Developed and made in Germany

magnetic ^emotion

Туре	luMIX	
Stirring points	1	
Stirring volume/stirring point	1 - 3,000 ml	
Speed range	100 - 2,000 rpm	
Stirring power (max.)	20 W	
Power setting	25/50/75/100% (4-step)	
Light setting	25/50/75/100% (4-step)	
Material housing	stainless steel	
Permitted operation conditions	-10 °C up to +40 °C (at 80% humidity)	
Electrical data	100-240 V / 50-60 Hz / 1.5 A	
Dimensions (WxDxH)	188 x 238 x 44 mm	
Protection category	IP65	
Weight (gross)	approx. 4.4 kg	
Order no.	31001	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Stirring - 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept for powerful mixing of solutions, extremely wide speed range from 100 to 2,000 rpm, jerk-free stirring even at low speeds, 4-step power setting for high power with regard to large and viscous stirring volumes and reduced power for the warming-free continuous operation.

Light - Bright and warming-free luminescent PMMA plate, white LEDs with homogenous brightness and variable light intensity (4-step), non-reflecting illumination..

Clear digital display for light-, speed- and power settings, SoftStart procedure for reliable catching/centering and safe accelerating of the stirring bar, QuickSet for quick setting of start and maximum speed, automatic memory function to save the last settings, outstanding long-life cycle of the hermetically sealed drive system, resistant stainless steel housing for easy and fast cleaning.

Accessories

ASTEROID 40 - Stirring bar Stirring bar remover PTFE

2mag MIXphilosophy

MIXdrive + MIXcontrol = unbeaten resistant and efficient stirring system



		ing systems	2mag - MIXphilosophy Combination possibilities of stirring systems					
Stirring drive	+	control unit						
MIXdrive 60 MIXdrive 15 MIXdrive 6 MIXdrive 1 MIXdrive 1 eco	MIXcontrol 40 (RS232) MIXcontrol 20 (RS232) (DINrail) MIXcontrol 20 (RS232) MIXcontrol eco (DINra							
cuvetteMIXdrive MIXdrive 60 HT								
MIXdrive 15 HT MIXdrive 6 HT	MIXcontrol 40 (RS232)	MIXcontrol 20 (RS232)	MIXcontrol eco (DINrail)					
MIXdrive 1 HT MIXdrive 1 eco HT	MIXcontrol 20 (RS232) MIXcontrol eco (DINrai							
	MIXdrive 15 MIXdrive 6 MIXdrive 1 MIXdrive 1 eco MIXdrive 1 XS cuvetteMIXdrive MIXdrive 60 HT MIXdrive 15 HT MIXdrive 6 HT MIXdrive 1 HT	MIXdrive 15 MIXdrive 15 MIXdrive 6 MIXdrive 1 MIXdrive 1 eco MIXdrive 1 eco MIXdrive 1 XS cuvetteMIXdrive MIXdrive 60 HT MIXdrive 60 HT MIXdrive 61 HT MIXdrive 1 S HT MIXdrive 1 S HT MIXdrive 1 S HT MIXdrive 1 S HT MIXdrive 1 HT MIXdrive 1 HT MIXdrive 1 eco HT	MIXdrive 15 MIXdrive 15 MIXdrive 6 MIXdrive 1 MIXdrive 1 MIXdrive 1 MIXdrive 1 MIXdrive 1 MIXdrive 1 MIXdrive 1 MIXdrive 1 MIXdrive 6 MIXdrive 6 MIXdrive 6 MIXdrive 6 MIXdrive 15 HT MIXcontrol 40 (RS232) MIXdrive 15 HT MIXcontrol 40 (RS232) MIXdrive 1 S MIXcontrol 20 (RS232) MIXdrive 1 S MIXcontrol 20 (RS232) MIXdrive 1 S MIXcontrol 20 (RS232) MIXdrive 1 S MIXcontrol 20 (RS232) MIXcontrol 20 (RS232)					

magnetic ^emotion

The individual combination possibilities of three control unit types MIXcontrol with the stirring drives MIXdrive always offer you the maximum quality for the optimum value for money exactly complementary to your needs.

The control units MIXcontrol define thereby the performance of the stirrers and the operation comfort.

The stirring drives MIXdrive are the indestructible equipment for your tasks in all days' laboratory work. No matter whether you decide for a serial device or a custom-made device, our aim is always: quality, stability and a reliable and affordable solution for exactly your needs. All MIXdrive are, with regard to the inner life, technically identical and can be combines with any of the 3 different control units MIXcontrol.

We offer you different stirring combinations for your individual applications. In case there will be no adequate stirrer available for you, we will be able to shortly work out a solution according to your needs.

Ask us – test us – we are there for you!



STIRRING DRIVES - TEMPERATURE RESISTANT AND SUBMERSIBLE

Designed for chemical, biotechnical and medical applications of all kinds, developed for maximum requirements, maximum mechanical, thermal and chemical stability as well as uncompromising continuous operation.



MIXdrive 1

Ultra flat, wear-free magnetic stirrer with 1 stirring position, suitable for vessels up to 10,000 ml.

- ℵ Wear-free magnetic stirrer, up to +200 °C
- ≈ Submersible, up to +95 °C
- ℰ Integrable in systems
- ℵ Developed and made in Germany





Туре	MIXdrive 1	
Stirring points	1	
Stirring volume/stirring point	1 - 10,000 ml	
Stirring power (max.)	20 W	
Material housing	stainless steel	
Perm. operation conditions (Basic)	-10 °C up to +50 °C in air, up to +50 °C in water	
Perm. operation conditions (HT)	up to +200 °C in air, up to +95 °C in water	
Operating voltage (max.)	48 VDC	
Dimensions (WxDxH)	180 x 180 x 38 mm	
Protection category	IP68	
Weight (gross)	approx. 3.5 kg	
Order no. Basic-version	40001	
Order no. HT-version	40201	

Individual units and dimensions on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

1 stirring position, 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept for powerful mixing of solutions, jerk-free stirring even at low speeds.

Fully encapsulated and hermetically sealed stainless steel housing, IP68, water-, dust-, and germproof, extremely robust construction and performance, easy to clean and maximum resistance, submersible in water- or oil baths, suitable for the use in CO₂-incubators and ovens.

Two different temperature areas available: **Basic-version** up to +50 °C, **HT-version** up to +95 °C submerged in water baths and up to +200 °C in ovens, individual dimensions on request.

Please order control unit separately!

Compatible control units

MIXcontrol eco MIXcontrol eco DINrail MIXcontrol 20 MIXcontrol 40 (only in combination with distriBOX 4)

Accessories

ASTEROID 40 - Stirring bar Stirring bar remover PTFE Extension cord for MIXdrive distriBOX 4 - distributor

Designed for chemical, biotechnical and medical applications of all kinds, developed for maximum requirements, maximum mechanical, thermal and chemical stability as well as uncompromising continuous operation.



MIXdrive 1 eco

Ultra flat, wear-free magnetic stirrer with 1 stirring position, suitable for vessels up to 3,000 ml.



- ≈ Submersible, up to +95 °C

- ℵ Developed and made in Germany



		Ama	Type Stirrin Stirrin Mater Perm. Opera Dimer Protec Weigh Order Order
			We reser

Туре	MIXdrive 1 eco	
Stirring points	1	
Stirring volume/stirring point	1 - 3,000 ml	
Stirring power (max.)	10 W	
Material housing	stainless steel	
Perm. operation conditions (Basic)	-10 °C up to +50 °C in air, up to +50 °C in water	
Perm. operation conditions (HT)	up to +200 °C in air, up to +95 °C in water	
Operating voltage (max.)	48 VDC	
Dimensions (WxDxH)	120 x 120 x 35 mm	
Protection category	IP68	
Weight (gross)	approx. 1.6 kg	
Order no. Basic-version	40101	
Order no. HT-version	40301	

Individual units and dimensions on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

1 stirring position, 100% maintenance-free and wear-free by inductive 2mag Magnetic-Drive concept for powerful mixing of solutions, jerk-free stirring even at low speeds.

Fully encapsulated and hermetically sealed stainless steel housing, IP68, water-, dust-, and germproof, extremely robust construction and performance, easy to clean and maximum resistance, submersible in water- or oil baths, suitable for the use in CO₂-incubators and ovens.

Two different temperature areas available: **Basic-version** up to +50 °C, **HT-version** up to +95 °C submerged in water baths and up to +200 °C in ovens, individual dimensions on request.

Please order control unit separately!

Compatible control units

MIXcontrol eco MIXcontrol eco DINrail MIXcontrol 20 (only in combination with distriBOX 4) MIXcontrol 40 (only in combination with distriBOX 4/8)

Accessories

ASTEROID 25 - Stirring bar Stirring bar remover PTFE Extension cord for MIXdrive distriBOX 4/8 - distributor

Designed for chemical, biotechnical and medical applications of all kinds, developed for maximum requirements, maximum mechanical, thermal and chemical stability as well as uncompromising continuous operation.



MIXdrive 1 XS

Ultra flat, wear-free magnetic stirrer with 1 stirring position, optimized for stirring volume up to 1,000 ml, ideal size for integration.

- ⋧ Wear-free magnetic stirrer, up to +200 °C
- ≈ Submersible, up to +95 °C
- ֎ Extremely robust construction
- ֎ Integrable in systems
- Neveloped and made in Germany





Туре	MIXdrive 1 XS		
Stirring points	1		
Stirring volume/stirring point	1 - 1,000 ml		
Stirring power (max.)	10 W		
Material housing	stainless steel		
Perm. operation conditions (Basic)	-10 °C up to +50 °C in air, up to +50 °C in wate		
Perm. operation conditions (HT)	up to +200 °C in air, up to +95 °C in water		
Operating voltage (max.)	48 VDC		
Dimensions (WxDxH)	48 x 48 x 18 mm		
Protection category	IP68		
Weight (gross)	approx. 0.36 kg		
Order no. Basic-version	40300		
Order no. HT-version	40303		

Individual units and dimensions on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

1 stirring position, 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept for powerful mixing of solutions, jerk-free stirring even at low speeds.

Fully encapsulated and hermetically sealed stainless steel housing, thread M4 integrated in stainless-steel bottom plate (not HT-version), IP68, water-, dust- and germproof, extremely robust construction and performance, easy to clean and maximum resistance, submersible in water- or oil baths, suitable for the use in CO_2 -incubators and ovens.

Two different temperature areas available: **Basic-version** up to +50 °C, **HT-version** up to +95 °C submerged in water baths and up to +200 °C in ovens, individual dimensions on request.

Please order control unit separately!

Compatible control units

MIXcontrol eco MIXcontrol eco DINrail MIXcontrol 20 (only in combination with distriBOX 4)

Accessories

ASTEROID 25 - Stirring bar Stirring bar remover PTFE Extension cord for MIXdrive distriBOX 4/8 - distributor

Designed for chemical, biotechnical and medical applications of all kinds, developed for maximum requirements, maximum mechanical, thermal and chemical stability as well as uncompromising continuous operation.



cuvetteMIXdrive 1

Miniaturized, wear-free magnetic stirrer with 1 stirring position for cuvettes, optimized for stirring volume up to 25 ml, ideal size for system integration e.g. in photometer.

- ℵ Maintenance- and wear-free
- ⋧ Submersible up to +50 °C
- ∂ Ideal for integration
- ∂ Outstanding long-life quality
- ֎ Developed and made in Germany

magnetic ^emotion



Туре	cuvetteMIXdrive 1	
Stirring points	1	
Stirring volume/stirring point	1 - 25 ml	
Stirring power (max.)	10 W	
Material housing	PUR	
Permitted operation conditions	-10 °C up to +50 °C in air, up to +50 °C in water	
Operating voltage (max.)	24 VDC	
Dimensions (WxDxH)	12.5 x 12. x 6 mm	
Protection category	IP68	
Weight (gross)	approx. 0.14 kg	
Order no.	40400	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

1 stirring position, 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept for reliable mixing of solutions, jerk-free stirring even at low speeds.

Fully encapsulated and hermetically sealed polyurethane housing, IP68, water-, dust- and germproof.

Extremely robust construction and performance, easy to clean and maximum resistance, submersible in water- or oil baths, suitable for the use in CO_2 -incubators and ovens.

Please order control unit separately!

Compatible control units

MIXcontrol eco MIXcontrol eco DINrail

Accessories

Extension cord for MIXdrive

Designed for reliable mass screenings as well as chemical, biotechnical and medical applications of all kinds, developed for the hard daily lab use and uncompromising continuous operation.



MIX 6 / MIX 15

Ultra-flat magnetic stirrer with 6 stirring positions, suitable for 6 x 1,000 ml Erlenmeyer flasks, or with 15 stirring positions, suitable for 15 x 250 ml beaker glasses (tall form).

- *№* Maintenance- and wear-free
- № For different vessel sizes
- ∂ Outstanding long-life quality
- ∂ Developed and made in Germany





Туре	MIX 6	MIX 15	
Stirring points	6	15	
Stirring point distance	130 mm	65 mm	
Stirring volume/stirring point	1 - 3,000 ml		
Speed range	100 - 2,	000 rpm	
Stirring power (max.)	20 W		
Power setting	5/10/15/20 W (4-step)		
Material housing	stainless steel		
Permitted operation conditions	-10 °C up to +50 °C (at 80% humidity)		
Electrical data	100-240 V / 50-60 Hz / 1.5 A		
Dimensions (WxDxH)	245 x 425 x 38 mm		
Protection category	IP65		
Weight (gross)	approx. 9.0 kg		
Order no.	30006 30015		

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

6 or 15 stirring positions, 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept, power reserves up to 3,000 ml stirring volume per stirring point, extremely wide speed range of 100 to 2,000 rpm, 100% synchronized speed, jerk-free stirring even at low speeds, 4-step power setting of the stirrer, high power setting for viscous liquids and reduced stirring power for long-term operation without any heating effects caused by the stirrer.

Clear digital display for settings of stirrer speed and power, SoftStart procedure for reliable catching/centering and safe acceleration of the stirring bar, QuickSet for quick setting of start- and maximum speed, Automatic memory function to save the last settings, outstanding long-life cycle of the hermetically sealed drive system, resistant stainless steel housing for easy and fast cleaning...

Accessories

ASTEROID 40 - Stirring bar Stirring bar remover PTFE

Designed for chemical, biotechnical and medical applications of all kinds, developed for the daily lab use and continuous operation.



MIX 15 eco

Ultra flat, wear-free magnetic stirrer with 15 stirring positions, suitable for 15 x 250 ml beaker glasses (tall form).

- ℵ Price-optimized, maintenance- and wear-free
- ℵ For different vessel sizes
- ֎ Developed and made in Germany

magnetic ^emotion



Price-optimized magnetic stirrer with 15 stirring positions, 100% maintenance-free and wear-free by inductive 2mag Magnetic-Drive concept for powerful mixing of solutions, power reserves up to 1,500 ml per stirring position, speed range from 120 up to 1,200 rpm, 100% synchronized speeds, jerk-free stirring even at low speeds. Clear digital display for stirrer speed settings, SoftStart procedure for reliable catching/centering and safe acceleration of the stirring bar. Automatic memory function to save the last settings, outstanding long-life cycle of the hermetically sealed drive system, resistant stainless steel housing for easy and fast cleaning.

Accessories

ASTEROID 40 - Stirring bar Stirring bar remover PTFE

Designed for reliable mass screenings as well as chemical, biotechnical and medical applications of all kinds, developed for the hard daily lab use and uncompromising continuous operation.



MIX 8 XL / MIX 12 XL

Ultra-flat magnetic stirrer, stirring point distance 90 mm, with 8 stirring positions, suitable for 8 x 600 ml beaker glasses (tall form), or with 12 stirring positions, suitable for 12 x 600 ml beaker glasses (tall form).

- ℵ New developed magnetic stirrer
- ℵ Maintenance- and wear-free
- ➢ For different vessel sizes
- ∂ Outstanding long-life quality
- 浸 Developed and made in Germany





Туре	MIX 8 XL	MIX 12 XL	
Stirring points	8	12	
Stirring point distance	90	mm	
Stirring volume/stirring point	1 - 3,000 ml		
Speed range	100 - 1,600 rpm		
Stirring power (max.)	20 W		
Power setting	5/10/15/20 W (4-step)		
Material housing	stainless steel		
Permitted operation conditions	-10 °C up to +50 °C (at 80% humidity)		
Electrical data	100-240 V / 50-60 Hz / 1.5 A		
Dimensions (WxDxH)	185 x 420 x 38 mm	275 x 420 x 38 mm	
Protection category	IP65		
Weight (gross)	approx. 7.8 kg	approx. 10.2 kg	
Order no.	30208	30212	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Powerful 8 or 12 stirring positions, 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept, power reserves up to 3,000 ml per stirring point, extremely wide speed range of 100 to 1,600 rpm, 100% synchronized speed, jerk-free stirring even at low speeds, 4-step power setting of the stirrer, high stirring power for viscous liquids and reduced stirring power for long-term operation without any heating effects caused by the stirrer.

Clear digital display for settings of stirrer speed and power, SoftStart procedure for reliable catching/centering and safe acceleration of the stirring bar, QuickSet for quick setting of start- and maximum speed, automatic memory function to save the last settings, outstanding long-life cycle of the hermetically sealed drive system, resistant stainless steel housing for easy and fast cleaning.

Accessories

ASTEROID 25 - Stirring bar Stirring bar remover PTFE

Designed for reliable mass screenings as well as chemical, biotechnical and medical applications of all kinds, developed for the hard daily lab use and uncompromising continuous operation.



MIX 4 MS

Ultra-flat magnetic stirrer with 4 individually controlled stirring positions, suitable for 4 x 2,000 ml Erlenmeyer flasks.

- ∂ 4 independent stirring positions
- ∂ Wear- and maintenance-free
- ⋧ Space-saving
- ∂ Outstanding long-life quality
- ∂ Developed and made in Germany

magnetic ^emotion



уре	MIX 4 MS	
tirring points	4	
tirring point distance	140 mm	
tirring volume/stirring point	1 - 3,000 ml	
peed range	100 - 2,000 rpm	
tirring power (max.)	28 W (4 x 7 W)	
ower setting	-	
1aterial housing	stainless steel	
ermitted operation conditions	-10 °C up to +50 °C (at 80% humidity)	
lectrical data	100-240 V / 50-60 Hz / 1.5 A	
imensions (WxDxH)	245 x 425 x 38 mm	
rotection category	IP65	
Veight (gross)	approx. 9.1 kg	
Irder no.	30404	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Powerful 4 stirring positions, 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept, power reserves up to 3,000 ml per stirring point, individual and independent speed setting for each stirring drive (Master-Slave-Technology), also ON/OFF function, synchronous speed mode for all positions, extremely wide speed range of 100 to 2,000 rpm, jerk-free stirring even at low speeds. Clear digital display for settings of stirrer speeds, SoftStart procedure for reliable catching/centering and safe a cceleration of the stirring bar, automatic memory function to save the last settings, outstanding long-life cycle of the hermetically sealed drive system, resistant stainless steel housing for easy and fast cleaning.

Accessories

ASTEROID 40 - Stirring bar Stirring bar remover PTFE

2mag MIXphilosophy

MIXdrive + MIXcontrol = unbeaten resistant and efficient stirring system



2mag - MIXphilosophy Combination possibilities of stirring systems				
	Stirring drive	+	control unit	
	MIXdrive 60			
	MIXdrive 15	MIXcontrol 40 (RS232)	MIXcontrol 20 (RS232)	MIXcontrol eco (DINrail)
Basic	MIXdrive 6	(=)	((2
up to +50 °C (in air)	MIXdrive 1	MIXcontrol	20 (RS232)	MIXcontrol eco (DINrail)
up to +50 °C (submersed)	MIXdrive 1 eco	MIXcontrol eco (DINrail)		
	MIXdrive 1 XS			DINrail)
	cuvetteMIXdrive			
	MIXdrive 60 HT			
НТ	MIXdrive 15 HT	MIXcontrol 40 (RS232)	MIXcontrol 20 (RS232)	MIXcontrol eco (DINrail)
(High Temperature)	MIXdrive 6 HT			
up to +200 °C (in air)	MIXdrive 1 HT	MIXcontrol	20 (RS232)	MIXcontrol eco (DINrail)
up to +95 °C (submersed)	MIXdrive 1 eco HT	MIXcontrol eco (DINrail)		
	MIXdrive 1 XS HT			

magnetic ^emotion

The individual combination possibilities of three control unit types MIXcontrol with the stirring drives MIXdrive always offer you the maximum quality for the optimum value for money exactly complementary to your needs.

The control units MIXcontrol define thereby the performance of the stirrers and the operation comfort.

The stirring drives MIXdrive are the indestructible equipment for your tasks in all days' laboratory work. No matter whether you decide for a serial device or a custom-made device, our aim is always: quality, stability and a reliable and affordable solution for exactly your needs. All MIXdrive are, with regard to the inner life, technically identical and can be combines with any of the 3 different control units MIXcontrol.

We offer you different stirring combinations for your individual applications. In case there will be no adequate stirrer available for you, we will be able to shortly work out a solution according to your needs.

Ask us – test us – we are there for you!



STIRRING DRIVES - TEMPERATURE RESISTANT AND SUBMERSIBLE

MULTIPLE STIRRING DRIVES WITH EXTERNAL CONTROL UNIT

Designed for chemical, biotechnical and medical applications of all kinds, developed for maximum requirements, maximum mechanical, thermal and chemical stability as well as uncompromising continuous operation.



MIXdrive 6 / MIXdrive 15 / MIXdrive 60

Ultra-flat, 100% maintenance- and wear-free magnetic stirrer with inductive 2mag-Magnetic-Drive concept for powerful mixing of solutions.

- ≈ Wear-free magnetic stirrer, up to +200 °C
- ≈ Submersible, up to +95 °C
- *№* Extremely robust construction
- *№* Integrable in systems
- Ś Developed and made in Germany





Туре	MIXdrive 6	MIXdrive 15	MIXdrive 60
Stirring points	6	15	60
Stirring point distance	130 mm	65 mm	32.5 mm
Stirring volume/stirring point	1 - 3.0	000 ml	1 - 500 ml
Stirring power (max.)	40 W		
Material housing	stainless steel		
Perm. operation conditions (Basic)	-10 °C up to +50 °C in air, up to +50 °C in water		
Perm. operation conditions (HT)	up to +200 °C in air, up to +95 °C in water		
Operating voltage (max.)	48 VDC		
Dimensions (WxDxH)	245 x 375 x 38 mm		
Protection category	IP68		
Weight (gross) approx.	9.4 kg	9.5 kg	8.9 kg
Order no. Basic-version	40006	40015	40060
Order no. HT-version	40206	40215	40260

Individual units and dimensions on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

MIXdrive 6

6 stirring positions, suitable for 6 x 1,000 ml Erlenmeyer flasks, stirring point distance 130 mm.

MIXdrive 15

15 stirring positions, suitable for 15 x 250 ml beaker glasses (tall form), stirring point distance 65 mm.

MIXdrive 60

60 (6x10) stirring positions, suitable for 60 testing tubes, ideal for integration, stirring point distance 32,5 mm.

100% synchronized speeds, jerk-free stirring even at low speeds. Fully encapsulated and hermetically sealed stainless steel housing, IP68, water-, dust-, and germproof.

Extremely robust construction and performance, easy to clean and maximum resistance, submersible in water- or oil baths, suitable for the use in CO₂-incubators and ovens.

Two different temperature areas available: Basic-version up to +50 °C, HT-version up to +95 °C submerged in water baths and up to +200 °C in ovens, individual dimensions on request.

Please order control unit separately!

Compatible control units

MIXcontrol eco MIXcontrol eco DINrail MIXcontrol 20 MIXcontrol 40

Accessories

ASTEROID 25/40 - Stirring bar Stirring bar remover PTFE Extension cord for MIXdrive



MIX control eco

Price-optimized control unit for small medium volumes and the connection of 1 stirring drive MIXdrive.



ℵ Speed range 120 up to 1,200 rpm

ℵ SoftStart

magnetic ^emotion

400 600 200 1000 off 1200	
man MIXcontrol eco	
1	

Туре	MIXcontrol eco		
Speed range	120 - 1,200 rpm		
Stirring power (max.)	10 W		
Power setting	-		
Material housing	plastic		
Perm. operation conditions	0 °C up to +40 °C (at 80% humidity)		
Electrical data	100-240 V / 50-60 Hz / 0.75 A		
Dimensions (WxDxH)	65 x 50 x 45 mm		
Protection category	IP20		
Weight (gross)	approx. 0.5 kg		
Order no.	90100		

Individual units on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Control unit with 10 watts stirring power, speed range from 100 up to 1,200 rpm, SoftStart for reliable catching/centering and safe accelerating of the stirring bar.

Compatible stirring drives

MIXdrive 1 MIXdrive 1 eco MIXdrive 1 XS cuvetteMIXdrive 1 MIXdrive 6 MIXdrive 15 MIXdrive 60



MIXcontrol eco DINrail

Price-optimized control unit for DIN installation (electrical cabinets).

- ➢ DINrail installation (electrical cabinets)
- 冷 Speed range 120 up to 1,200 rpm
- NoftStart





Туре	MIXcontrol eco DINrail		
Speed range	120 - 1,200 rpm		
Stirring power (max.)	10 W		
Power setting	-		
Material housing	plastic		
Perm. operation conditions	0 °C up to +40 °C (at 80% humidity)		
Operating voltage (max.)	24 VDC		
Dimensions (WxDxH)	24 x 100 x 120 mm		
Protection category	IP20		
Weight (gross)	approx. 0.25 kg		
Order no. potentiometer	90140		
Order no. 0-10 V	90150		
Order no. RS232	90160		

Individual units on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Control unit for DINrail installation (electrical cabinets) for small to medium volumes and the connection of one stirring drive MIXdrive, with analog interface 0-10 V, RS232 or potentiometer for ON/OFF and speed adjustment. Control unit with 10 watts stirring power, speed range from 100 to 1,200 rpm, SoftStart for reliable catching/centering and safe accelerating of the stirring bar.

Compatible stirring drives

MIXdrive 1 MIXdrive 1 eco MIXdrive 1 XS cuvetteMIXdrive 1 MIXdrive 6 MIXdrive 15 MIXdrive 60



MIXcontrol 20

Universal control unit for the connection of 1 stirring drive MIXdrive.

- ≈ 20 watts stirring power
- ⋧ 4-step power setting
- ℵ Speed range 100 1,600 rpm
- ֎ Developed and made in Germany

magnetic ^emotion





Туре	MIXcontrol 20	
Speed range	100 - 1,600 rpm	
Stirring power (max.)	20 W	
Power setting	25/50/75/100% (4-step)	
Material housing	stainless steel	
Permitted operation conditions	0 °C up to +40 °C (at 80% humidity)	
Electrical data	100-240 V / 50-60 Hz / 1.5 A	
Dimensions (WxDxH)	200 x 155 x 38 mm	
Protection category	IP20	
Weight (gross)	approx. 1.9 kg	
Order no.	90200	
Order no. RS232	90210	

Individual units on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Control unit with 20 watts stirring power, speed range from 100 up to 1,600 rpm, jerk-free stirring also at low speeds, power setting (4-steps) for high power with regard to large and viscous stirring amounts and reduced power for the warming-free continuous operation of MIXdrive stirring drives.

Clear digital display for speed- and power settings, SoftStart for reliable catching/centering and safe accelerating of the stirring bar. QuickSet for quick setting of start- and maximum speed, automatic memory function to save the last settings, resistant stainless steel housing, easy to clean.

Also available with RS232 interface.

Interfaces (0-10 V, 4-20 mA, USB, Ethernet, OPC UA, LADS oriented), other housings e.g. for DINrail installation and customized software adaption on request.

Compatible stirring drives

MIXdrive 1

MIXdrive 1 eco (only in combination with distriBOX 4) MIXdrive 1 XS (only in combination with distriBOX 4) MIXdrive 6 MIXdrive 15 MIXdrive 60



MIX control 40

Powerful control unit, for 2mag MIXdrive series, developed for the hard daily lab use and uncompromising continuous operation.

- ℵ 40 watts stirring power
- ➢ For connection of 2 stirring drives MIXdrive
- ⋧ Speed range 100 2,000 rpm
- ֎ Developed and made in Germany







Туре	MIXcontrol 40	
Speed range	100 - 2,000 rpm	
Stirring power (max.)	40 W	
Power setting	10 - 100% (10-step)	
Material housing	stainless steel	
Permitted operation conditions	0 °C up to +40 °C (at 80% humidity)	
Electrical data	100-240 V / 50-60 Hz / 1.5 A	
Dimensions (WxDxH)	200 x 155 x 38 mm	
Protection category	IP20	
Weight (gross)	approx. 1.9 kg	
Order no.	90400	
Order no. RS232	90410	

Individual units on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Powerful control unit for the connection of up to two stirring drives MIXdrive, powerful 40 watts stirring power, wide speed range from 100 up to 2,000 rpm, jerk-free stirring also at low speeds, power setting (10-steps) for high power with regard to large and viscous stirring amounts and reduced power for the warming-free continuous operation of MIXdrive stirring drives. Clear digital display for speed- and power settings, SoftStart for reliable catching/centering and safe accelerating of the stirring bar. QuickSet for quick setting of start- and maximum speed, automatic memory function to save the last settings, resistant stainless steel housing, easy to clean.

Also available with RS232 interface.

Interfaces (0-10 V, 4-20 mA, USB, Ethernet, OPC UA, LADS oriented), other housings e.g. for DINrail installation and customized software adaption on request.

Compatible stirring drives

MIXdrive 1 (only in combination with distriBOX 4) MIXdrive 1 eco (only in combination with distriBOX 4/8) MIXdrive 1 XS (only in combination with distriBOX 4/8) MIXdrive 6 MIXdrive 15 MIXdrive 60

HEATABLE STIRRER

Developed for stirring and temperating jobs in daily lab use. Designed for chemical, biotechnical and medical applications of all kinds, developed for the hard daily lab use and uncompromising continuous operation.



hotMIX 1

Smallest hotplate magnetic stirrer of the world. Compact, powerful and very robust magnetic stirrer with aluminium hotplate, unbeaten space-saving and also suitable for the operation with viscous and aggressive media.

- ⋧ Fast electrical heating
- ℵ Robust, long-life quality
- ֎ Developed and made in Germany

magnetic ^emotion

Туре	hotMIX 1
Stirring points	1
Stirring volume/stirring point	1 - 2,000 ml
Speed range	100 - 1,600 rpm
Stirring power (max.)	10 W
Power setting	-
Temperature range	+40 °C up to +200 °C
Material heating plate	aluminium alloy
Material housing	stainless steel
Permitted operation conditions	-10 °C up to +40 °C (at 80% humidity)
Operating voltage	230 V
Heating power (max.) at 230 V	500 W
Dimensions (WxDxH)	120 x 140 x 80 mm
Protection category	IP20
Weight (gross)	approx. 2.0 kg
Order no.	61001

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Price-optimized magnetic stirrer, 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept for mixing of solutions, speed range from 100 up to 1,600 rpm, jerk-free stirring even at low speeds. SoftStart for reliable catching/centering and safe accelerating of the stirring bar, resistant stainless steel housing for easy and fast cleaning, spill over protection, extremely robust construction. Hotplate made from resistant aluminium alloy, max. heating temperature +200 °C, socket for contact thermometer (accessory), warning-LED in case of hot plate, 3 integrated overtemperature protection loops against overheating.

Accessories

mixTHERM - Thermometer

For mass screenings, incubating or concentrating of probes, evaporating of solvents, simple and easy system integration.



STIRRING HOTPLATE 6 / STIRRING HOTPLATE 15

Stirring-heating plate with 6 stirring positions, stirring point distance 130 mm, suitable for 6 x 1,000 ml Erlenmeyer flasks or with 15 stirring positions, stirring point distance 65 mm, suitable for 15 x 250 ml beaker glasses (tall form).

- ⋧ Fast heating-up
- ≈ Energy-saving
- ℵ For all vessel sizes and shapes
- ⋧ Simple system integration
- Developed and made in Germany





Туре	HOTPLATE 6	HOTPLATE 15	
Stirring points	6 15		
Stirring point distance	130 mm 65 mm		
Stirring volume/stirring point	1 - 1,500 ml	1 - 1,000 ml	
Stirring power (max.)	40 W		
Power setting	10 - 100% (10-step)		
Temperature range	+40 °C up to +200 °C		
Speed range	100 - 2,000 rpm		
Operating voltage	230 V		
Heating power (max.) at 230 V	1,200 W		
Dimensions (WxDxH)	250 x 430 x 75 mm		
Protection category	IP20		
Weight (gross)	approx. 13 kg		
Order no.	66306	66315	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Stirring - Powerful 6 or 15 stirring points, maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept, extremely wide speed range of 100 to 2,000 rpm, 100% synchronized speed, jerk-free stirring even at low speeds, 10-step power setting of the stirrer, high power setting for viscous liquids and reduced stirring power for long-term operation without any heating effects caused by the stirrer, clear digital display for settings of stirrer speed, stirrer power and temperature of the integrated heater, SoftStart procedure for reliable catching/centering and safe acceleration of the stirring bar. **Heating** - Heating plate made of resistant aluminium alloy, PTFE-coated for increased chemical resistance and easier cleaning, fast heating-up by powerful integrated electrical heater, homogenous temperature distribution, maximum heating temperature +200 °C, PowerDetect for automatic compensation of physically determined loss in stirring power which can come up to 40% at high temperatures, independent over-temperature protection inside the plate. Inclusive separate control unit for operation of magnetic stirrer and integrated heater, for long-term protection of the electronics against aggressive liquids and gases as well as thermal influences, stainless steel housing.

Accessories

External probe sensor Pt100 Stand for heatMIXcontrol

STIRRING-HEATING BLOCKS WITH EXTERNAL CONTROL UNIT

For viscosity measurements, R&D, combinatorial chemistry, biological applications, testing the solution behavior of synthetic materials, quality control in food, chemical and waste water technology, simple and easy system integration.



STIRRING DRYBATH 15-100 / STIRRING DRYBATH 15-250

Stirring-heating block combination with 15 bore holes, suitable for 100 ml lab-flasks and for 150 ml beaker glasses tall form (ø 57.5 mm, depth 48 mm, distance 65 mm) or with 15 bore holes suitable for 250 ml lab-flasks and for 400 ml beaker glasses tall form (ø 71.5 mm, depth 78 mm, distance 84 mm).

- ∂ Fast heating-up
- ֎ Insulated frame
- Norable quality
- Developed and made in Germany

Ty S1 S1 S1 S1 S1 S1 S1 C Te S1 O H D P1 W O





уре	DRYBATH 15-100	DRYBATH 15-250	
tirring points	15		
tirring point distance	65 mm	84 mm	
Bore holes Ø/T	Ø 57.5 / 48 mm	Ø 71.5 / 78 mm	
tirring volume/stirring point	1 - 150 ml	1 - 400 ml	
tirring power (max.)	40 W		
Power setting	10 - 100% (10-step)		
emperature range	+40 °C up to +200 °C		
peed range	100 - 2,000 rpm		
Dperating voltage	230 V		
leating power (max.) at 230 V	1,000 W	1,150 W	
Dimensions (WxDxH)	280 x 430 x 120 mm	335 x 515 x 150 mm	
Protection category	IP20		
Veight (gross)	approx. 20 kg	approx. 35 kg	
Order no.	63100	63250	

Individual units and dimensions on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Stirring - maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept, extremely wide speed range of 100 to 2,000 rpm, 100% synchronized speed, jerk-free stirring even at low speeds, 10-step power setting of the stirrer, high stirring power for viscous liquids and reduced stirring power for long-term operation without any heating effects caused by the stirrer, clear digital display for settings of stirrer speed, stirrer power and temperature of the integrated heater, SoftStart procedure for reliable catching/ centering and safe acceleration of the stirring bar. **Heating** - Heating plate made of resistant aluminium alloy, PTFE-coated for increased chemical resistance and easier cleaning, fast heating-up by powerful integrated electrical heater, homogenous temperature distribution, maximum heating temperature +200 °C, PowerDetect for automatic compensation of physically determined loss in stirring power which can come up to 40% at high temperatures, independent over-temperature protection inside the plate. Inclusive separate control unit for operation of magnetic stirrer and integrated heater, for long-term protection of the electronics against aggressive liquids and gases as well as thermal influences, stainless steel housing.

Accessories

External probe sensor Pt100 Stand for heatMIXcontrol Reducing adapters - Inserts for reducing the bore hole diameter

STIRRING-HEATING BLOCKS WITH EXTERNAL CONTROL UNIT

For viscosity measurements, R&D, combinatorial chemistry, biological applications, testing the solution behavior of synthetic materials, quality control in food, chemical and waste water technology, simple and easy system integration.



STIRRING DRYBATH 8-250 ERL

Stirring-heating block combination with 8 stirring positions (ø 87.2 mm, depth 78 mm, distance 100/140 mm), suitable for 250 ml Erlenmeyer flasks.

- ⋧ Fast heating-up
- ℵ Insulated frame
- ⋧ Durable quality
- ֎ Developed and made in Germany

Type



DRYBATH 8-250 ERL



Stirring points	8		
Stirring point distance	100 / 140 mm		
Bore holes Ø/T	Ø 87.2 / 78 mm		
Stirring volume/stirring point	1 - 250 ml		
Stirring power (max.)	40 W		
Power setting	10 - 100% (10-step)		
Temperature range	+40 °C up to +200 °C		
Speed range	100 - 2,000 rpm		
Operating voltage	230 V		
Heating power (max.) at 230 V	1,150 W		
Dimensions (WxDxH)	335 x 515 x 150 mm		
Protection category	IP20		
Weight (gross)	approx. 40 kg		
Order no.	63825		

Individual units and dimensions on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Stirring - Maintenance-free and wear-free by inductive 2mag-Magnetic Drive concept, extremely wide speed range of 100 to 2,000 rpm, 100% synchronized speed, jerk-free stirring even at low speeds, 10-step power setting of the stirrer, high stirring power for viscous liquids and reduced stirring power for long-term operation without any heating effects caused by the stirrer, clear digital display for settings of stirrer speed, stirrer power and temperature of the integrated heater, SoftStart procedure for reliable catching/centering and safe acceleration of the stirring bar. **Heating** - Massive heating block made of resistant aluminium alloy, PTFE-coated for increased chemical resistance and easier cleaning, contact adapters for an optimal heat transfer, lowest possible temperature gradient inside the stirring vessels, integrated electrical heater, maximum heating temperature +200 °C.

PowerDetect for automatic compensation of physically determined loss in stirring power which can come up to 40% at high temperatures, specially made insulated stainless steel frame for energy- and time saving heating process, additionally for a homogenous temperature distribution and protection against accidental contact, independent over-temperature protection inside the plate. Inclusive separate control unit for operation of magnetic stirrer and integrated heater, for long-term protection of the electronics against aggressive liquids and gases as well as thermal influences, stainless steel housing.

Inclusive 8 contact adapters for an optimized heat transfer.

Accessories

External probe sensor Pt100 Stand for heatMIXcontrol

STIRRER FOR CELL CULTURES WITH INTERNAL CONTROL

Designed for careful and protective mixing of cell suspensions and culture broths, developed for highest demands as well as warming-free and uncompromising continuous operations.



bioMIX 1

Price-optimized maintenance-free magnetic stirrer with 1 stirring point, large surface, space saving, also suitable for viscous culture media.

- ⋧ Space saving
- ℵ Germ-proof, easy to clean
- Outstanding long-life quality
- ֎ Developed and made in Germany

magnetic ^emotion



Туре	bioMIX 1	
Stirring points	1	
Stirring volume/stirring point	5 - 5,000 ml	
Speed range	5 - 250 rpm	
Stirring power (max.)	6 W	
Power setting	10 - 100% (10-step)	
Material housing	stainless steel	
Permitted operation conditions	-10 °C up to +50 °C (at 80% humidity)	
Electrical data	100-240 V / 50-60 Hz / 1.5 A	
Dimensions (WxDxH)	188 x 238 x 38 mm	
Protection category	IP64	
Weight (gross)	approx. 3.3 kg	
Order no.	85001	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Maintenance-free 2mag-stepper-motor-drive for extremely slow and 100% jerk-free and gentle mixing of cell cultures, extended speed range from 5 up to 250 rpm, jerk-free, even and smooth stirring also at slow speeds, power setting (10-step) for high power with regard to large and viscous stirring amounts and reduced power for the warming-free continuous operation. Digital display for speed- and power setting, SoftStart for reliable and safe accelerating of stirring bar, QuickSet for quick setting of start- and maximum speed, saving of last working status when turning off, suitable for culture flasks with spinning propeller- and spinning ball systems.

Fully encapsulated and water resistant stainless steel housing, germ-proof, IP64, easy to clean and firm, robust construction and performance, ultra-flat design.

Hint

For using in incubators we recommend due to the minimized temperature emission the especially developed series bioMIXdrive.

STIRRING DRIVES FOR CELL CULTURES WITH EXTERNAL CONTROL UNIT

Designed for careful and protective mixing of cell suspensions and culture broths, developed for highest demands as well as warming-free and uncompromising continuous operations for e.g. in CO₂-incubators.



bioMIXdrive 1 / 2 / 3 / 4

Ultra flat, wear-free magnetic stirrer with large standing surface, up to 5,000 ml stirring volume, also suitable for viscous culture media.



- ⋧ Warming-free long term operation in CO₂-incubator
- ℵ Outstanding long-life quality
- ℵ Developed and made in Germany

Type



bioMIXdrive



11				
Stirring points	1	2	3	4
Stirring point distance	- 140 mm			
Stirring volume/stirring point	5 - 5,000 ml			
Speed range	5 - 250 rpm			
Stirring power/stirring point		5	W	
Material housing	stainless steel			
Perm. operation conditions	-10 °C up to +50 °C (at 100% humidity)			
Perm. operation conditions	+50 °C submerged in water			
Operating voltage (max.)	48 VDC			
Dimensions (WxD, mm)	180 x 180	130 x 270	130 x 410	270 x 270
Dimensions (H, mm)	38			
Protection category	IP68			
Weight (gross), approx.	2.5 kg	2.9 kg	4.1 kg	5.6 kg
Order no.	80001	80002	80003	80004

Individual units and dimensions on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Maintenance-free 2mag-stepper-motor-drive for extremely slow and 100% jerk-free and gentle mixing of cell cultures, extended speed range from 5 up to 250 rpm, jerk-free, even and smooth stirring also at slow speeds, suitable for culture flasks with spinning propeller- and spinning ball systems.

Fully encapsulated and hermetically closed stainless steel housing, IP68, water-, dust- and germ-proof, extremely robust construction and performance, easy to clean and maximum firm. Submersible in water baths, designed for the warming-free long term operation in CO_2 -incubators, individual measurements on request.

Please order control unit separately!

Compatible control units

bioMIXcontrol bioMIXcontrol S bioMIXcontrol 4MS (only for bioMIXdrive 1)

Accessories

Extension cord for bioMIXdrive

CONTROL UNITS FOR bioMIXdrive STIRRING DRIVES

Designed for bioMIXdrive stirring drives. For chemical, biotechnical and medical applications of all kinds.



bioMIXcontrol / bioMIXcontrol S

Efficient control unit, for 2mag bioMIXdrive series, developed for the hard daily lab use and uncompromising warming-free continuous operation. Also available as stackable control unit bioMIXcontrol S for system integration.

- ℵ Speed range 5 250 rpm
- ℵ Power setting 10-step
- ℵ Outstanding long-life quality
- ℵ Developed and made in Germany

magnetic ^emotion





Туре	bioMIXcontrol	bioMIXcontrol S	
Speed range	5 - 250 rpm		
Stirring power (max.)	40 W		
Power setting	10 - 100% (10-step)		
Material housing	stainless steel		
Permitted operation conditions	0 °C up to +40 °C (at 80% humidity)		
Electrical data	100-240 V / 50-60 Hz / 1.5 A		
Dimensions (WxDxH)	200 x 155 x 38 mm 215 x 120 x 57 mm		
Protection category	IP20		
Weight (gross)	approx. 1.9 kg		
Order no.	98400	98500	

Individual units on request. 3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Control unit for the connection of 1 stirring drive bioMIXdrive 1, 2, 3 or 4 extended speed range from 5 to 250 rpm, without any own-generated heating effect, jerk-free stirring, even and smooth stirring also at slow speeds.

Power setting (10-steps) for high power with regard to large and viscous stirring amounts and reduced power for the warming-free continuous operation e.g. in CO_2 -incubators.

Digital display for speed- and power setting, SoftStart for reliable and safe accelerating of stirring bar, QuickSet for quick setting of start- and maximum speed, storage of last working status when turning off, firm stainless steel housing, easy to clean.

Interfaces (0-10 V, 4-20 mA, RS232, USB, Ethernet, OPC UA, LADS oriented), other housings e.g. for DINrail installation and firmware adaptions on request.

Compatible stirring drives

bioMIXdrive 1 bioMIXdrive 2 bioMIXdrive 3 bioMIXdrive 4



bioMIXcontrol 4MS

Efficient control unit, for 4 drives 2mag bioMIXdrive 1, developed for the hard daily lab use and uncompromising warming-free continuous operation.

- ℵ Speed range 5 250 rpm
- ➢ For 4 stirring drives bioMIXdrive 1
- ∂ Outstanding long-life quality
- ℵ Developed und Made in Germany

magnetic ^emotion



We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Control unit for the connection of 4 stirring drive bioMIXdrive 1, individual and independent speed setting for each stirring drive (Master-Slave-Technology), also ON/OFF function, synchronous speed mode for all drives, extended speed range from 5 to 250 rpm, 6 watts/stirring position without any own-generated heating effect, automatic stirring power for the warming-free continuous operation e.g. in CO₂-incubators, jerk-free stirring, even and smooth stirring also at slow speeds. Digital display for speed setting, SoftStart for reliable and safe accelerating of stirring bar, QuickSet for quick setting of start and maximum speed, storage of last working status when turning off, firm stainless steel housing, easy to clean. Compatible stirring drives bioMIXdrive 1

MICROTITER PLATE STIRRERS

Designed for reliable mass screenings as well as chemical, biotechnical and medical applications of all kinds, developed for the hard daily lab use and uncompromising continuous operation.



MIXdrive 6 / 12 / 24 / 96 MTP

Ultra flat, wear-free magnetic stirrer with 6/12/24/96 stirring positions, suitable for microtiter plates.

- *№* Wear-free magnetic stirrer
- ⋧ Submersible, up to + 50 °C, integrable
- ➢ Extremely robust construction
- ֎ Developed und Made in Germany





Type MIXdrive	6 MTP	12 MTP	24 MTP	96 MTP
Stirring points	6	12	24	96
Stirring point distance	39 mm	26 mm	19 mm	9 mm
Stirring volume/stirring point (max.)	20 ml	7 ml	3,5 ml	0,5 ml
Stirring power	10 W			
Material housing	stainless steel			
Permitted operation conditions	-10 up to +50 °C in air, up to +50 °C in water submerged			
Operating voltage	48 VDC			
Dimensions (WxDxH)	140 x 102 x 26 mm			
Protection category	IP68			
Weight (gross)	approx. 1.3 kg			
Order no.	43006	on request	43024	43096

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

All stirring positions 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept for powerful mixing of solutions, 100% synchronized speeds, jerk-free stirring even at low speeds.

Fully encapsulated and hermetically sealed stainless steel housing, IP68, water-, dust- and germproof, extremely robust construction and performance, easy to clean and maximum resistance, submersible in water- resp. oil baths up to +50°C. Suitable for the use in CO_2 incubators and ovens, removable centering pins, suitable for integration into robots.

Please order control unit separately.

Compatible control units

MIXcontrol eco MIXcontrol MTP

CONTROL UNITS FOR MIXDRIVE MTP STIRRING DRIVES

Designed for MIXdrive MTP stirring drives.



MIX control MTP

Powerful control unit for connection of one stirrring drive MIXdrive 6/12/24/96 MTP.

- ➢ For stirring drives MIXdrive MTP
- ⋧ Extremely robust construction
- ℵ Developed und Made in Germany

magnetic ^emotion



Universal control unit for the connection of one stirring drive MIXdrive MTP, universal 10 watts stirring power, wide speed range from 100 up to 2,000 rpm, jerk-free stirring also at low speeds, 4-step power setting.

Clear digital display for speed- and power settings, SoftStart for reliable catching/centering and safe accelerating of the stirring bar.

QuickSet for quick setting of start- and maximum speed, automatic memory function to save the last settings, resistant stainless steel housing, easy to clean.

Also available with RS232 interface.

Interfaces (0-10 V, 4-20 mA, USB, Ethernet, OPC UA, LADS oriented), other housings e.g. for DINrail installation and customized software adaption on request.

Compatible stirring drives

MIXdrive 6 MTP MIXdrive 12 MTP MIXdrive 24 MTP MIXdrive 96 MTP

INDUSTRIAL STIRRERS FOR HIGHLY VISCOUS MEDIA AND LARGE VOLUMES

Designed for chemical and biotechnical applications of all kinds as e.g. within the field of quality insurance and pilot plant stations, developed for highest demands, unbeaten strong, high stability as well as uncompromising continuous operation.



MIX 1 XL

Ultra flat, strong magnetic stirrer with large set-up area and fully automatic stirring bar monitoring mixWATCH, up to 40 litres stirring volume, also suitable for highly viscous media, very strong magnetic coupling, therefore predestined for magnetic power transmission over wide distances or for thick vessel bottoms.

- ∂ Up to 40 litres or for viscous media
- *∂* mixWATCH stirring bar monitoring
- ∂ Outstanding long-life quality
- ∂ Developed and made in Germany

magnetic ^emotion



Туре	MIX 1 XL	
Stirring points	1	
Stirring volume/stirring point	1 - 40 litres	
Speed range	70 - 2,000 rpm	
Stirring power (max.)	20 W	
Power setting	5/10/15/20 W (4-step)	
Acceleration	variable, 20/30/60/120 s	
Interval	variable, 0 - 240 s (2-step)	
Material housing	stainless steel	
Permitted operation conditions	-10 °C up to +50 °C (at 95% humidity)	
Electrical data	100-240 V / 50-60 Hz / 1.5 A	
Dimensions (WxDxH)	225 x 280 x 38 mm	
Protection category	IP64	
Weight (gross)	approx. 4.5 kg	
Order no.	30201	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Maintenance-free 2mag-motor-drive with large speed range from 70 up to 2,000 rpm, fully encapsulated and tight stainless steel housing (IP64), extremely robust construction and performance, easy to clean and firm.

Integrated controller with strong 20 watts stirring power, 4-step power setting for high power with regard to large and viscous stirring volumes and reduced power for the warming-free continuous operation, SoftStart-function (adjustable, 4 steps) for reliable catching/centering and safe accelerating of the stirring bar at different volumes and viscosities, adjustable interval times for automated restart of the stirrer after max. 240 seconds.

Digital display for speed-, power-, SoftStart- and interval setting, QuickSet for quick setting of start and maximum speed, after-run and glass breakage, storage of last function status when turning off.

mixWATCH - Stirring bar monitoring with fully automatic trim of the maximum reachable speed for a reliable and unattended operation of the stirrer (detailed description on page 39).

Accessories

ASTEROID 40 - Stirring bar SATELLITE 70 - Beared sirring bar Stirring bar remover PTFE

INDUSTRIAL STIRRERS FOR HIGHLY VISCOUS MEDIA AND LARGE VOLUMES

Designed for chemical and biotechnical applications of all kinds as e.g. within the field of quality insurance and pilot plant stations, developed for highest demands, unbeaten strong, high stability as well as uncompromising continuous operation.



max MI X

Unbeaten strong magnetic stirrer with large set-up area, up to 150 litres stirring volume, also suitable for highly viscous media.



- ℵ Compact dimensions
- ℵ Ideal for highly viscous media
- ֎ Developed and made in Germany

magnetic ^emotion

Туре
Stirring
Stirring
Speed
Stirring
Power
Acceler
Materia
Permitt
Electric
Dimens
Protect
Weight
Order r
We reserve

2	maxMIX	
ing points	1	
ing volume/stirring point	1 - 150 litres	
ed range	100 - 1,200 rpm	
ing power (max.)	40 W	
er setting	-	
eleration	variable, 20 - 99 s	
erial housing	stainless steel	
nitted operation conditions	-10 °C up to +50 °C (at 80% humidity)	
trical data	100-240 V / 50-60 Hz / 4.2 A	
ensions (WxDxH)	265 x 320 x 68 mm	
ection category	IP64	
ght (gross)	approx. 10.0 kg	
er no.	30501	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Very strong magnetic coupling, therefore predestined for magnetic power transmission over wide distances (up to 60 mm), for thick vessel bottoms or for double-coat vessels. Maintenance-free 2mag-motor-drive with large speed range from 100 up to 1,200 rpm, fully encapsulated and tight stainless steel housing, IP64, extremely robust construction and performance, easy to clean under running water and firm. Integrated control unit with strong 40 watts stirring power, large speed range from 100 up to 1,200 rpm, SoftStart-function (variable 20 - 99 seconds) for reliable catching/centring and safe accelerating of the stirring bar at different volumes and viscosities.

Digital display for speed- and SoftStart-setting, QuickSet for quick setting of the start and maximum speed, QuickStopfunction against after-run and glass breakage, storage of last function status when turning off, firm stainless steel housing, easy to clean.

Accessories

ASTEROID 70 - Stirring bar SATELLITE 140 - Beared stirring bar for large volumes Stirring bar remover XL maxFRAME - Enlargement of surface area

INDUSTRIAL STIRRING SYSTEMS FOR HIGHLY VISCOUS MEDIA AND LARGE VOLUMES

Designed for chemical and biotechnical applications of all kinds as e.g. within the field of quality insurance and pilot plant stations, developed for highest demands, unbeaten strong, high stability as well as uncompromising continuous operation.



MIXdrive 1 XL

Ultra flat, compact and strong magnetic stirrer, up to 40 litres stirring volume, also suitable for highly viscous media, temperature resistant up to +100°C.

MIXcontrol XL

Powerful control unit for MIXdrive 1 XL stirring drives with 20 watts stirring power and mixWATCH stirring bar monitoring.

- ℵ Temperature resistant up to 100°C
- ℵ Up to 40 litres stirring volume
- ⋧ mixWATCH stirring bar monitoring
- ֎ Outstanding long-life quality
- Neveloped and made in Germany



Туре	MIXdrive 1 XL	MIXcontrol XL
Stirring points	1	-
Stirring volume/stirring point	1 - 40 litres	-
Stirring poweer (max.)	-	20 W
Power setting	-	5/10/15/20 W (4-step)
Speed range	-	70 - 2,000 rpm
Acceleration	-	variable, 20/30/60/120 s
Interval	-	variable, 0 - 240 s
Material housing	stainless steel	
Operating voltage (max.)	24 VDC	-
Electrical data	-	100-240 V / 50-60 Hz / 1.5 A
Permitted operation conditions	+5 °C up to +100°C (at 95% humidity)	0 °C up to +40°C (at 80% humidity)
Dimensions (WxDxH)	225 x 225 x 38 mm	200 x 155 x 38 mm
Protection category	IP66	IP20
Weight (gross)	approx. 4.0 kg	approx. 1.9 kg
Order no.	40601	90600

Individual units and dimensions on request.

3 years warranty on material and manufacturing.

We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Maintenance-free 2mag-motor-drive with large speed range from 70 up to 2,000 rpm, IP66, temperature resistant up to +100°C, designed for use in ovens, extremely robust construction and performance, easy to clean and firm, very strong magnetic coupling by SmCo-magnets, therefore predestined for magnetic power transmission over wide distances or for thick vessel bottoms.

The stirring system MIXdrive 1 XL with the control unit MIXcontrol XL is equipped as standard with **mixWATCH**. (detailed description on page 39).

Please order control unit separately!

SoftStart-function (adjustable, 4-step), for reliable catching/ centering and safe accelerating of the stirring bar at different volumes and viscosities, adjustable interval times for automated restart of the stirrer after max. 240 seconds.

20 watts stirring power, speed range from 70 up to 2,000 rpm, power setting in 4 steps. Digital display for speed and interval setting, QuickSet for quick setting of start and maximum speed, QuickStop-function against after-run and glass breakage, storage of last function status when turning off. Equipped as standard with RS232 interface.

Accessories

ASTEROID 40 - Stirring bar SATELLITE 70 - Beared sirring bar Stirring bar remover PTFE Extension cord for MIXdrive 1 XL

Interfaces (0-10 V, 4-20 mA, USB, Ethernet, OPC UA, LADS oriented), other housings e.g. for DINrail installation and customized software adaption on request.
INDUSTRIAL STIRRING SYSTEMS FOR HIGHLY VISCOUS MEDIA AND LARGE VOLUMES

Designed for chemical and biotechnical applications of all kinds as e.g. within the field of quality insurance and pilot plant stations, developed for highest demands, unbeaten strong, high stability as well as uncompromising continuous operation.



MAXdrive

Unbeaten strong magnetic stirrer with large set-up area, up to 250 litres stirring volume, also suitable for highly viscous media.

MAXcontrol

Control unit for the connection of MAXdrive, powerful 60 watts stirring power , large speed range from 100 up to 1,200 rpm.

- ℵ Up to 250 litres stirring volume
- ✤ Stirring over wide distances
- ℵ Ideal for highly viscous media
- Outstanding long-life quality
- Developed and made in Germany





Туре	MAXdrive MAXcontrol		
Stirring points	1	-	
Stirring volume/stirring point	1 - 250 litres	-	
Stirring power (max.)	-	60 W	
Power setting	-	-	
Speed range	-	100 - 1,200 rpm	
Acceleration	-	variable, 20 - 99 s	
Material housing	stainless steel		
Operating voltage (max.)	24 VDC	-	
Electrical data	-	100-240 V / 50-60 Hz / 4.2 A	
Permitted operation conditions	-10 °C up to +50°C (at 95% humidity)	0 °C up to +40°C (at 80% humidity)	
Dimensions (WxDxH)	260 x 260 x 68 mm	200 x 175 x 51 mm	
Protection category	IP64	IP20	
Weight (gross)	approx. 9.8 kg	approx. 2.0 kg	
Order no.	50001	94400	

Individual units and dimensions on request. 3 years warranty on material and manufacturing.

We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Very strong magnetic coupling, therefore predestined for magnetic power transmission over wide distances (up to 60 mm), for thick vessel bottoms or for double-coat vessels, ideal for the integration into process engineering systems.

Maintenance-free 2mag-motor-drive with large speed range from 100 up to 1,200 rpm. Fully encapsulated and tight stainless steel housing, IP64, extremely robust construction and performance, easy to clean under running water and maximum firm. Individual measurements on request.

Please order control unit separately!

Variable SoftStart-function for reliable catching/centering and safe accelerating of the stirring bar at different volumes and viscosities. Digital display for speed- and SoftStart-setting (variable 20 - 99 seconds).

QuickSet for quick setting of the start-and maximum speed, QuickStop-function against after-run and glass breakage, storage of last function status when turning off, firm stainless steel housing, easy to clean.

Interfaces (0-10 V, 4-20 mA, RS232, USB, Ethernet, OPC UA, LADS oriented), other housings e.g. for DINrail installation and customized software adaption on request.

Accessories

ASTEROID 70 - Stirring bar SATELLITE 140 - Beared stirring bar for large volumes Stirring bar remover XL Extension cord for MAXdrive/FABdrive maxFRAME - Enlargement of surface area

INDUSTRIAL STIRRING SYSTEMS FOR HIGHLY VISCOUS MEDIA AND LARGE VOLUMES

Designed for chemical and biotechnical applications of all kinds as e.g. within the field of quality insurance and pilot plant stations, developed for highest demands, unbeaten strong, high stability as well as uncompromising continuous operation.



FABdrive

Unbeaten strong magnetic stirrer with very compact housing dimensions, up to 1,000 litres stirring volume, also suitable for highly viscous media.

FABcontrol

Control unit for the connection of magnetic stirrer FABdrive, powerful 100 watts stirring power, large speed range from 80 up to 990 rpm.

- ℵ Up to 1,000 litres stirring volume
- ℰ Stirring over wide distances
- ֎ Ideal for highly viscous media
- ℵ Developed and made in Germany



We

Туре	FABdrive	FABcontrol	
Stirring volume/stirring point	1 - 1,000 litres	-	
Speed range	-	80 - 990 rpm	
Stirring power (max.)	-	100 W	
Power setting	-	-	
Acceleration	-	variable, 20 - 99 s	
Buzzer, Tolerance	-	variable, 5 - 50 rpm	
Material housing	stainless steel		
Perm. operation conditions	-10 °C up to +50 °C (at 95% humidity)	0 °C up to +40 °C (at 80% humidity)	
Operating voltage (max.)	24 VDC	-	
Electrical data	-	100-240 V / 50-60 Hz / 4.2 A	
Dimensions (WxDxH)	180 x 180 x 75 mm	200 x 175 x 51 mm	
Protection category	IP64	IP20	
Weight (gross)	approx. 4.8 kg	approx. 2.0 kg	
Order no.	50101	94402	

Individual units and dimensions on request.

3 years warranty on material and manufacturing. Ve reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing

Very strong magnetic coupling, therefore predestined for magnetic power transmission over wide distances (up to 150 mm), for thick vessel bottoms or for double-coat vessels, ideal for the integration into process engineering systems and industrial processing.

Maintenance-free 2mag-motor-drive with large speed range from 80 up to 990 rpm. Fully encapsulated and tight stainless steel housing, IP64, extremely robust construction and performance, easy to clean under running water and maximum firm. Individual measurements on request.

Please order control unit separately!

Variable SoftStart-function for reliable catching/centering and safe accelerating of the stirring bar at different volumes and viscosities. Digital display for speed- and SoftStart-setting (variable, 20 - 99 seconds) and live speed, QuickSet for quick setting of the start-and maximum speed, QuickStop-function against after-run and glass breakage.

Beep-function at falling below set speed, storage of last function status when turning off, firm stainless steel housing, easy to clean.

Accessories

ASTEROID 70 - Stirring bar SATELLITE 140 - Beared stirring bar for large volumes Stirring bar remover XL Extension cord for MAXdrive/FABdrive fabFRAME - Enlargement of surface area

Interfaces (0-10 V, 4-20 mA, RS232, USB, Ethernet, OPC UA, LADS oriented), other housings e.g. for DINrail installation and customized software adaption on request.



2mag mixWATCH

Stirring bar monitoring with fully automatic trim of the maximum reachable speed for a reliable and unattended operation of the stirrer.

≈ Contactless

Network Automatic

➢ Integrated in MIX 1 XL, MIXdrive 1 XL, steriMIXdrive





An absolutely innovation is the integrated contact-free stirring bar control mixWATCH. In case of the magnetic coupling is separated between the magnetic stirrer and the stirring bar during the acceleration phase e.g. caused by to high set speed or changing viscosities, the magnetic stirrer will be stopped, the stirring bar centred and started again. For the magnetic stirres MIX 1 XL and MIXdrive 1 XL, the set speed level will be reduced automatically by 100 rpm to ensure a reliable stirring process.

The stirring drive steriMIXdrive starts also automatically, but without reducing the stirring speed.

mixWATCH is ensuring a very high process safety, e.g. with using of closed non-glass vessels or at unattended overnight operation.

SPECIAL STIRRING SYSTEM FOR OPERATION IN PRESSURE, VACUUM AND STEAM

Designed for operation in pressure, vacuum and steam up to +150°C, e.g. in steam sterilizers.



steriMIXdrive

Ultra flat, compact and strong magnetic stirrer, up to 40 litres stirring volume, also suitable for use in autoclaves.

steriMIXcontrol

Powerful control unit with 20 watts stirring power and mixWATCH stirring bar monitoring (without stirring speed reduction).

Very strong magnetic coupling by SmCo-magnets, therefore predestined for magnetic power transmission over wide distances or for thick vessel bottoms. Maintenance-free 2mag-motor-drive with large speed range from 100 up to 1,200 rpm.

Fully encapsulated, pressure- and vacuum resistant stainless steel housing, IP68, temperature resistant up to +150°C, extremely robust construction and performance, easy to clean and firm.

easyMOUNT plug coupling for quick installation of the stirrer cable through a pressure and steam tight grommet into a steam sterilizer.

Please order control unit separately!

SoftStart-function, for reliable catching/centering and safe accelerating of the stirring bar at different volumes and viscosities, adjustable interval times for automated restart of the stirrer after max. 240 sec.

Digital display for speed- and interval setting, QuickSet for quick setting of start and maximum speed, QuickStopfunction against after-run and glass breakage, storage of last function status when turning off.

- ℵ Up to 40 litres stirring volume
- ➢ mixWATCH stirring bar monitoring
- ℰ Steam-, pressure- and vakuum resistant
- Ideal für high viscous media
- Neveloped and made in Germany



Туре	steriMIXdrive	steriMIXcontrol
Stirring points	1	-
Stirring volume/stirring point	1 - 40 litres	-
Stirring power (max.)	-	20 W
Power setting	-	-
Speed range	-	100 - 1,200 rpm
Acceleration	-	variable, 20/30/60/120 s
Interval	-	variable, 0 - 240 s
Material housing	stainless steel / PPSU hermetically closed	stainless steel
Perm.operation conditions	-10 °C up to +150 °C (at 100% humidity)	0 °C up to +40 °C (at 80% humidity)
Operating voltage (max.)	24 VDC	-
Electrical data	-	100-240 V / 50-60 Hz / 1.5 A
Dimensions (WxDxH)	140 x 140 x 32 mm	200 x 155 x 38 mm
Protection category	IP68	IP20
Weight (gross)	approx. 1.5 kg	approx. 1.9 kg
Order no.	40700	95200

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Accessories

ASTEROID 40 - Stirring bar SATELLITE 70 - Beared sirring bar Stirring bar remover PTFE

Interfaces (0-10 V, 4-20 mA, RS232, USB, Ethernet, OPC UA, LADS oriented), other housings e.g. for DINrail installation and customized software adaption on request.

The stirring system steriMIXdrive with the control unit steriMIXcontrol is equipped as standard with **mixWATCH**. (detailed description on page 39).

STIRRING DRIVE FOR EXPLOSION PROTECTION ZONES

Designed for chemical, biotechnical and medical applications of all kinds, developed for the daily lab use and continuous operation in explosion protection zones.



atexMIXdrive

Explosion protected, ultra flat, wear-free magnetic stirrer with 1 stirring position, for up to 10,000 ml.

- ⋧ Explosion protected magnetic stirring system
- ℵ Wear-free magnetic stirrer
- ≈ Up to 10,000 ml strirring volume
- ℵ Outstanding long-life quality
- ֎ Developed and made in Germany





Туре	atexMIXdrive
Stirring points	1
Stirring volume/stirring point	1 - 10,000 ml
Stirring power (max.)	20 W
Material housing	stainless steel
Permitted operation conditions	-10 °C up to +40 °C in air, up to +40 °C in water
Operating voltage (max.)	48 VDC
Dimensions (WxDxH)	180 x 180 x 38 mm
Protection category	IP68
Weight (gross)	3.5 kg
Ex-mark	ATEX II 2G Ex mb IIC T4 Gb
Order no.	49001

2 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

1 stirring position, 100% maintenance-free and wear-free by inductive 2mag-Magnetic-Drive concept for powerful mixing of solutions, jerk-free stirring even at low speeds.

Fully encapsulated and hermetically sealed stainless steel housing, IP68, water-, dust- and germproof. Extremely robust construction and performance, easy to clean and maximum resistance. Submersible in water- or oil baths, suitable for the use in CO_2 -incubators and ovens, temperate range up to +40 °C, certified according to ATEX II 2G Ex mb IIC T4 Gb.

Please order control unit separately!

Hint: Stirring drive atexMIXdrive and control unit atexMIXcontrol (cleanroom) are firmly connected!

Compatible control units

atexMIXcontrol Explosion protected control unit (protection zone 1) atexMIXcontrol cleanroom Explosion protected control unit (protection zone 2) MIXcontrol 20 Universal control unit (outside the protection zone)

Accessories

ASTEROID 40 - Stirring bar Stirring bar remover PTFE

CONTROL UNITS FOR atexMIXDRIVE STIRRING DRIVE

Designed for 2mag stirring drives atexMIXdrive.

Designed for chemical, biotechnical and medical applications of all kinds.



atexMIX control

Explosion protected control unit for the connection of 1 stirring drive atexMIXdrive, suitable for protection zone 1.

atexMIXcontrol cleanroom

Explosion protected control unit for the connection of 1 stirring drive atexMIXdrive, suitable for protection zone 2 or cleanrooms.

- ⋧ Explosion protected control unit
- ≈ 20 watts stirring power
- ≈ 4-step power setting
- ⋧ Speed range 100 2,000 rpm
- ֎ Developed and made in Germany





Туре	atexMIXcontrol	atexMIXcontrol cleanroom	
Speed range	100 - 2,00	00 rpm	
Stirring power (max.)	20 V	V	
Power setting	5/10/15/20	W (4-step)	
Material housing	aluminium/plastic	aluminium	
Perm. operation conditions	0 °C up to +40 °C (at 80% humidity)		
Electrical data	100-240 V / 50-60 Hz / 1.5 A		
Dimensions (WxDxH)	360 x 460 x 200 mm	210 x 55 x 30 mm	
Weight (gross)	approx. 18 kg	approx. 4.5 kg	
Protection category	IP65		
Ex-mark	ATEX II 2G Ex db eb IIC T6 Gb	ATEX II 3G Ex nA nL IIC T6	
Order no.	99200	99300	

2 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

20 watts stirring power, speed range from 100 up to 2,000 rpm, jerk-free stirring even at low speeds, 4-step power setting.

Clear digital display for speed- and power settings, SoftStart for reliable catching/centering and safe accelerating of the stirring bar.

QuickSet for quick setting of start and maximum speed, automatic memory function to save the last settings, explosion protected housing, operation temperature up to +40°C, certified according to ATEX II 2G Ex db eb IIC T6 Gb. 20 watts stirring power, speed range from 100 up to 2,000 rpm, jerk-free stirring even at low speeds, 4-step power setting.

Clear digital display for speed- and power settings, SoftStart for reliable catching/centering and safe accelerating of the stirring bar.

QuickSet for quick setting of start and maximum speed, automatic memory function to save the last settings, explosion protected housing, operation temperature up to +40°C, certified according to ATEX II 3G Ex nA nL IIC T6. Hint: Control unit atexMIXcontrol (cleanroom) and stirring drive atexMIXdrive are firmly connected!

Compatible stirring drives

atexMIXdrive Explosion protected stirring drive

STIRRING BARS

Designed for strong stirrers and viscous media. Designed for chemical, biotechnical and medical applications of all kinds.



ASTEROID 25 / 40 / 70

New developed, high-effective stirring bar, ideal for large mixing volumes, viscous media, stirring over wide distances or for disadvantageous vessel bottoms.





Туре	ASTEROID 25	ASTEROID 40	ASTEROID 70
Material (Cover)	PTFE		
Material (Magnet)	SmCo		
Shape	triangular, convex		
Dimensions (HxL)	14 x 25 mm 28 x 40 mm 39 x 70 mm		
Dimensions (Ø)	16 mm	33 mm	45 mm
Weight (gross) approx.	0.07 kg	0.12 kg	0.37 kg
Order no.	44025	44040	44070

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

More than 4x higher torque transmissible, unbeatable increased mixing effect in comparison to conventional stirring bars, high energetic and long-term stable magnetic field, no demagnetization effect, convex design for a minimum of bearing surface and reliable rotation, triangular shape for powerful shear force, better mixing effect at lower rotation speed. 300% less PTFE abrasion than comparable round shape stirring bars, significantly higher rotation speed reachable, high quality, even and sealed PTFE cover, FDA-conform, steam sterilizable at 121°C.

Multiple improved performances for all types of magnetic stirrers.

Accessories

Stirring bar remover PTFE Stirring bar remover XL

STIRRING BARS

Designed for strong stirrers and viscous media. Designed for chemical, biotechnical and medical applications of all kinds.



SATELLITE 70

New developed, high-effective stirring bar, recommended for magnetic stirrers MIXdrive 1 XL or MIX 1 XL.

SATELLITE 140

New developed, high-effective stirring bar, recommended for magentic stirrers FABdrive, MAXdrive or maxMIX.

magnetic ^emotion





Туре	SATELLITE 70	SATELLITE 140
Shape	round, tripod	round, tripod
Material (Cover)	PTFE	ZEDEX
Dimensions ØxL bars	Ø 13 x 70 mm	Ø 32 x 140 mm
Dimensions Ø tripod construction	Ø 99 mm	Ø 227 mm
Weight (gross)	approx. 0.15 kg	approx. 0.46 kg
Order no.	44700	44900

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Extremely strong, abrasion-free and wear-free magnetic stirring bar, bearing on a tripod-construction, ideal for large mixing volumes, viscous media, stirring over wide distances or for disadvantageous vessel bottoms, especially for the use in stainless steel vessels and in the case of media containing abrasive particles, e.g. sand or sediments.

No direct contact between stirring bar and reaction vessel, therefore steady firm, wear-free, stirring bar sealed with PTFE, tripod construction made of ZEDEX, chemical resistance comparable to PTFE, FDA-certification, wearing parts can be re-ordered separately. Extremely strong, abrasion-free and wear-free magnetic stirring bar, bearing on a tripod-construction, ideal for large mixing volumes, viscous media, stirring over wide distances or for disadvantageous vessel bottoms, especially for the use in stainless steel vessels and in the case of media containing abrasive particles, e.g. sand or sediments.

No direct contact between stirring bar and reaction vessel, therefore steady firm, wear-free, chemical resistance comparable to PTFE, FDA-certification, wearing parts can be re-ordered separately.

Accessories

Stirring bar remover PTFE Stirring bar remover XL

Designed for strong stirring bars.



Stirring bar remover PTFE

PTFE-coated flexible stirring bar remover with extremely strong permanent magnet (neodym), 70 cm length.

Stirring bar remover XL

New developed, extremely strong stirring bar remover with an integrated corrosion-resistant high energy magnet, ideal for reliable inserting and removing of strong and heavy magnetic stirring bars.



magnetic	emotion
----------	---------

Туре	Stirring bar remover PTFE	Stirring bar remover XL
Material	PTFE	stainless steel
Dimensions ØxL	Ø 15 x 700 mm	Ø 28 x 1,000 mm
Weight (gross)	approx. 0.5 kg	approx. 1.8 kg
Order no.	45070	45100

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Universal chemical resistance also in aggressive media, ideal for fast removal of heavy and big magnetic stirring bars with a weight up to 400 g.

Simple handling due to a special developed stripping mechanism, length of 100 cm for deep reaction vessels, construction of stainless steel, solid and easy to clean, high chemical resistance.

Ideal for fast removal of big and heavy magnetic stirring bars with a weight up to 1,000 g.

For 2mag stirring systems with external control.



Extension cords

Extension cord for stirring drives, length 3m.

distriBOX 4/8

4-fold / 8-fold distributor.

magnetic ^emotion

49008

MIXdrive MIXdrive 1 XL bioMIXdrive FABdrive/MAXdrive



Length			3 m	
Perm. operation condition	ns	-10°C up to +50°C (at 80% humidity)		
Weight (gross) approx.		0.15 kg		0.27 kg
Order no.	46100	46108	46200	46300
Туре	dist	riBOX 4		distriBOX 8
Connection sockets		4		8
Dimensions (WxDxH)	48 x 48	48 x 48 x 25 mm		x 92 x 25 mm
Weight (gross) approx.	0.	0.18 kg		0.27 kg

Extension cords for

Order no.

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

49004

Extension cord for MIXdrive, length 3m. Extension cord for MIXdrive 1 XL, length 3m. Extension cord for bioMIXdrive, length 3m. Extension cord for MAXdrive and FABdrive, length 3m. Distributor for synchronized operating of 4 or 8 stirring drives with one MIXcontrol 20 or MIXcontrol 40, identical operation conditions, synchronous speeds, same stirring power, space-saving, sealed stainless steel housing. Designed for 2mag multiple stirrers.



mixRACK

Stand for test tubes.

magnetic ^emotion



For exactly centred positioning of test tubes on the stirring points, rack function, can be used to carry all test tubes at once, two convenient handles, stainless steel construction. Designed for 2mag STIRRING DRYBATH and STIRRING HOTPLATE.



heatMIXcontrol - spare part

Separate control unit heatMIXcontrol for operation of STIRRING DRYBATH and STIRRING HOTPLATE.

Stand for heatMIXcontrol

Stand for vertical positioning of the control unit heatMIXcontrol, incl. assembling material.



magnetic ^emotion

Туре	heatMIXcontrol - spare part	
Stirring power (max.)	40 W	
Power setting	10 - 100% (10-step)	
Speed range	100 - 2,000 rpm	
Temperature range	+40 °C up to +200 °C	
Material housing	stainless steel	
Permitted operation conditions	0 °C up to +40 °C (at 80% humidity)	
Electrical data	230-240 V / 50-60 Hz / 5 A	
Dimensions (WxDxH)	225 x 171 x 48 mm	
Weight (gross)	approx. 2.2 kg	
Protection category	IP20	
Order no.	68300	
Туре	Stand heatMIXcontrol	
Material	stainless steel	
Dimensions (WxDxH)	225 x 132 x 150 mm	
Weight (gross)	approx. 1.0 kg	
Order no.	69500	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Long-term protection of the electronics against aggressive liquids and gases as well as thermal influences, independent and adjustable over-temperature limiter for the additional protection of the heating controller against technical defects or unintended operation errors, stainless steel housing, compact and space-saving, connection for external Pt100 sensor, RS232 interface. Stable stainless steel stand for space-saving positioning of the control unit heatMIXcontrol, organized cable route on rear side, easy installation.

Compatible stirring drives

STIRRING DRYBATH 15-100 STIRRING DRYBATH 15-250 STIRRING DRYBATH 8-250 ERL STIRRING HOTPLATE 6 STIRRING HOTPLATE 15 Designed for 2mag STIRRING DRYBATH and STIRRING HOTPLATE.



External probe sensor Pt100

External probe sensor Pt100 for control unit heatMIXcontrol.

Reducing adapter

Inserts (aluminium alloy) for STIRRING DRYBATH 15-100 and STIRRING DRYBATH 15-250 for reducing the bore hole diameter.

magnetic ^emotion



Туре	External probe sensor Pt100	
Sensor type	Pt100	
Material	stainless steel	
Dimensions (ØxL)	Ø 3 x 150 mm	
Cable length	1 m	
Weight (gross)	approx. 0.1 kg	
Order no.	69100	
Reducing adapter	for DRYBATH 15-100	for DRYBATH 15-250
Outer dimensions (ØxH)	Ø 57 5 x 47 mm	Ø 71 5 x 78 mm

Reducing adapter	for DRYBATH 15-100	for DRYBATH 15-250
Outer dimensions (ØxH)	Ø 57.5 x 47 mm	Ø 71.5 x 78 mm
Inner diameter	on request	
Order no.	on request	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Temperature sensor for direct controlling of the probe temperature.

Compatible control units heatMIXcontrol For reduction of the bore hole diameter, guarantees an excellent heat transfer in the case of using smaller glasses. Individual inner diameter on request.

Compatible stirring drives

STIRRING DRYBATH 15-100 Stirring-heating block combination, 15 stirring positions for 100 ml laboratory flasks.

STIRRING DRYBATH 15-250 Stirring-heating block, 15 stirring positions for 250 ml laboratory flasks. Designed for 2mag hotMIX.

Mag magnetic motion

mixTHERM

Thermometer with flexible Pt 1000 sensor, precise and fast rules by fuzzy logic control.

magnetic ^emotion



Туре	mixTHERM	
Measuring range	-50 °C up to +300 °C	
Weight (gross)	approx. 0.7 kg	
Order no.	61099	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Precise and fast control by Fuzzy-Logic-Technology, measuring range -50 °C up to +300 °C, accuracy +/- 1 °C, identification of probe breakdown and probe short circuit, safety- and control circuit according to DIN 12878 Class 1 and 2, display background illumination, LED indication for control status, IP 65, approx. 150 g, incl. clamp for support rod.

Compatible magnetic stirrer hotMIX Heatable stirrer, 1 stirring position

netic stirrer

Designed for large volumes.



maxFRAME

Frame for easy and fast enlargement of surface area and loading capacity of magnetic stirrers maxMIX and MAXdrive.

fabFRAME

Frame for easy and fast enlargement of surface area and loading capacity of magnetic stirrer FABdrive.

magnetic ^emotion



Туре	maxFRAME	
Material	stainless steel	
Dimensions (WxDxH)	500 x 500 x 70 mm	
Weight (gross)	approx. 10.5 kg	
Order no.	48500	
Туре	fabFRAME	
Material	stainless steel	
Dimensions (WxDxH)	500 x 500 x 76 mm	
Weight (gross)	approx. 10.5 kg	
Order no.	48600	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

Solid welded stainless steel construction, easy to clean due to enclosed design, high chemical resistance, simple enlargement of the utility space of maxMIX and MAXdrive to 500 x 500 mm, loading capacity up to 1000 kg.

No installation necessary.

Solid welded stainless steel construction, easy to clean due to enclosed design, high chemical resistance, simple enlargement of the utility space of FABdrive to 500 x 500 mm, loading capacity up to 1000 kg.

No installation necessary.



mixBATH

Durable, robust stainless steel bath, easy to clean, suitable fro MIXdrive 6/15/60 and MIXdrive 1/1 eco/1 XS.

mixBATHcover

Chemically resistant, transparent cover for mixBATH, consisting of polycarbonate, for minimization of evaporation effects.

magnetic ^emotion

40815

mixBATH



Material	stainless steel	
Iviaceriai	Stallliess Steel	
Inner dimensions (WxDxH)	277 x 537 x 175 mm	
Outer dimensions (WxDxH)	318 x 600 x 190 mm	
Weight (gross)	approx. 8.0 kg	
Order no.	40950	
mixBATHcover	for MIXdrive 6	for MIXdrive 15
Material	polycarbonate	
Dimensions (WxDxH)	315 x 575 x 8 mm	
Weight (gross)	approx. 2.6 kg	

3 years warranty on material and manufacturing. We reserve the right to make technical alterations. We do not assume liability for errors in printing or typeing.

40806

Fitting for immersion thermostats (heating/ cooling), easy to clean, even weld and electropolished, framed ridge formed as handle, integrated drain makes emptying liquid easy and safe, exact positioning of the stirrer in the case of use the stirrer under the bath. Stirring point mark inside the bath, left-/right stop for optimized thermostat positioning (circulation pump) for use with MIXdrive 6 or MIXdrive 15. Different cutouts for the perfect positioning of the thermostat in the case of using the stirring drives MIXdrive 6 or MIXdrive 15 immersed into the mixBATH.

Compatible stirring drives

MIXdrive 6 MIXdrive 15 MIXdrive 60

Type

Order no.

MIXdrive 1 MIXdrive 1 eco MIXdrive 1 XS



mixBATH PC

Durable, robust bath, made of polycarbonate, easy to clean, suitable for MIXdrive 6/15/60 and MIXdrive 1/1 eco/1 XS.





Туре	mixBATH PC	
Material	polycarbonate	
Inner dimensions (WxDxH)	280 x 540 x 178 mm	
Outer dimensions (WxDxH)	300 x 560 x 198 mm	
Weight (gross)	approx 8.3 kg	
Order no.	40850	

3 ans de garantie sur le matériau et la fabrication. Sous réserve de modifications techniques. Pas de responsabilité pour erreurs d'impression ou d'ortographe.

Water bath made of polycarbonate, durable up to +50° C, robust, easy to clean, suitable for MIXdrive 6/15/60 as well as and MIXdrive 1/1 eco/1 XS., fitting for immersion thermostats(heating/cooling), easy to clean, exact positioning of the stirrer in the case of use the stirrer under the bath.

Compatible stirring drives

MIXdrive 6 MIXdrive 15 MIXdrive 60

MIXdrive 1 MIXdrive 1 eco MIXdrive 1 XS

SPEZIALISED STIRRING SYSTEMS AND CUSTOMIZED STIRRES FOR INDIVIDUAL REQUIREMENTS



2mag - Customized Stirrers

The main importance of our passion lies upon our customers and our products!



magnetic ^emotion

Your desired stirrer - configuration options

- Individual housing dimensions
- Stirring positions from 1 to 96
- Individual arrangement of stirring positions
- Variable cable length and positioning of cable output
- Adapation to environmental conditions (temperature, pressure, steam, etc.)
- Control/Interfaces (RS232, 0-10 V, 4-20 mA)
- Other housings of control units e.g. for DINrail installation
- Special functions (Interval, Timer, etc.)
- Customized firmware (Default values)
- Integration (screwing, covering, assembly)
- Customized labeling (elektrolysis)

2mag is providing you more than 30 years of competence in the conception, development, production and sales of magnetic and inductive drive concepts, as e.g. magnetic stirrers.

This experience paired with a great passion and emotion for our customers – and our products – enables the fast and economic creation of solutions precisely for your needs. "Customizing", the manufacturing of special made products, is our business, our competence and our passion! Individual requests will result in individual – and even more important – also economically realizable solutions.

Ask us – test us – we are there for you!



2mag - Sustainability

The trend theme sustainability of 2mag is not only a short-living term of a current spirit of time but a concept and philosophy since the founding in 2007. At the beginning as we designed our products, we thought highly of a maximal reliability, quality, durability and low lifecycle costs.

magnetic ^emotion



Convince yourself of the sustainability and the value of 2mag magnetic stirrers!

- Longlife quality due to hermetically closed stainless steel housings, completely sealed
- 2mag magnetic stirring drives are water-, dust- and germ-proofed and even in some cases pressureand vacuum-tight
- High efficiency, 30 40% more stirring power than other comparable products of our competitors at the same electrical input

- 100% wear- and maintenance-free
- High quality Zero-stand-by power supply units for minimization of energy costs in the stand-by mode
- Regional suppliers warrant a high quality and a minimum environmental impact due to short transport routes
- Excellent price/performance with regard to trouble-free economic life of 10 years and more in daily use
- 3 years warranty





magnetic ^emotion 2mag AGITATION - the world is moving CERTIFIED I am interested in 2mag products. e-mail contact: 2mag AG Please contact me. Schragenhofstrasse 35 I-K phone contact: Company / Organisation 80992 Muenchen Germany Department Name / First name +49 (89) 38 15 31 10 Street info@2mag.de www.2mag.de Zip / Town Please e-mail to: info@2mag.de L Stamp 2mag is not only active and successful in the field of laboratory equipment. Also for the international famous light-design company Ingo Maurer in Munich we created a soundless magnetic drive for an exclusive illuminating and furnishing object. Further appliances not really typical for the laboratory field are to be developed by us with a lot of passion - be surprised!

2mag – your contact for sophisticated and also fulfillable solutions for your needs and issues in the range of mixing, tempering and dosing.

Delirium Yum

Light object, designed by Ingo Maurer and Sebastian Hepting www.ingo-maurer.com